

PRESTO W92x

Powerful Temperature Control Systems for demanding applications with a temperature range from -92 °C to +250 °C

PRESTO® W92 - Highly Dynamic Temperature Control Systems The PRESTO® W92 Highly Dynamic Temperature Control Systems are amongst the most powerful units available. Their impressive power is harnessed by cutting-edge control technology. The W92 provides extraordinary heating, cooling, and pumping performance. But it also comes with the functions and features that make the new generation of PRESTO® so unique.

Your advantages

- Powerful circulating pumps, electronically adjustable in stages or by setting the pressure value Choose between a magnetically coupled centrifugal pump or a magnetically coupled gear pump for pressures up to 5.5 bar and virtually constant flow rate at any pressure
- Integrated 5.7" industrial color touchscreen displays all essential information and enables simple fingertip control
- Extensive warning, protection, and monitoring functions with detailed self-explanatory messages
- ICC cascade control for extraordinary precision, temperature stability $\pm 0.05\text{ °C} \dots \pm 0.2\text{ °C}$
- Ambient temperature range $+5\text{ °C}$ to $+40\text{ °C}$
- Integrated programmer with real-time clock
- Filling level indicator and pump capacity displayed electronically
- Interface for SD memory card
- Connections for USB, Ethernet, RS232, and Alarm output
- Optional analog connections, RS485, Profibus DP, Modbus
- 2-stage water-cooled refrigeration system



Made in Germany

Technical Data

Order No.	9421923							
Category	Temperature Control PRESTO							
Working temperature range (°C)	-92 ... +250							
Temperature control	ICC							
Temperature stability (°C)	$\pm 0.05 \dots \pm 0.2$							
Setting / display resolution	0.01 °C							
Temperature Display	TFT Touchscreen							
Heating capacity (kW)	18							
Cooling capacity (Medium: JULABO Thermal Ethanol)	°C	200	20	0	-20	-40	-60	-80
	kW	29	25	18	11	10	7	1.5
Pump capacity flow rate (l/min)	18 ... 70							
Pump capacity flow pressure (psi)	11.6 ... 79.77							
Viscosity max. (cSt)	70							
Pump connections	M38x1.5							
Refrigerant	R23, R404A							

External Pt100 sensor connection	integrated
Digital interface	RS232, SD memory card, USB, Ethernet, Modbus, Alarm-out Optional: RS485, Profibus
Analog connection input / output	Optional
Ambient temperature	5 ... 40 °C
Dimensions W x L x H (inch)	37.4 x 50 x 74.8
Weight (LBS)	1764
Sound pressure level (distance 1 m) max. (dBA)	74
Process volume min. (active heat exchanger volume) liters	28 (16)
Internal usable expansion vol. (liters)	40
Classification according to DIN12876-1	Classification III (FL)
Cooling of compressor	2-stage Water
Cooling water connection	G ¾" male with barbed fittings for tubing ½" ID
Cooling water consumption (l/min)	16...43
Cooling water temperature (°C)	<30
Cooling water differential pressure (bar)	0.5
Power requirement V / Hz / A	3x 480/60/38
Available voltage versions	3 x 400V/50Hz (+/- 10%) Without Plug 3 x 480V/60Hz (+/- 10%) Without Plug

All data refers to the nominal voltage of 400 V, 3-phase, nominal frequency of 50 Hz and ambient temperature of +20 °C. Cooling capacity measured at max. pump stage. All pump data refers to a bath fluid with a specific density of 1 kg/dm³.

Tip: Counter-cooling your PRESTO with a Recirculating Cooler

If there is no cooling water, the PRESTO W92x can be cooled down with a recirculating cooler with a cooling capacity of 45 kW at a flow temperature of 15°C. The required circulating pump has to ensure a flow rate of 43 l/min at a counter-pressure of 0.5 bar. The recommended minimum tank volume is 200 liters.

Characteristics

Display

TFT DISPLAY **State-of-the-art display technology**
TFT Display for comfortable user guidance, colored display of measurement values, graphs and control options, user-defined views

Operation

TOUCH **Optimal ease of use**
Touch screen for direct operation via display

HELP **Instructions inside**
Help menus and explanations in plain text for all control options, help messages and warning messages

MULTI LINGUAL **Multilingual user guidance**
Language selection for display of control options, notifications and warning messages via touchscreen

ACCESS LEVELS **Convenience for several users**
Administrator level for customizing instrument settings, user levels with limited permissions for fast and safe defined access, password protection, all levels adjustable

Temperature Control

ICC **For perfect results**
'Intelligent Cascade Control', automatic & self optimizing adjustment of PID control parameters, temperature stability ±0.01 °C ... <±0.2 °C

TCF **Full control**
'Temperature Control Features', for individual optimization, access to all important control parameters, additional settings for band limit, limits, co-speedfactor etc.

Pt100 Control from the external application
 External Pt100 sensor connection for precise measurement and control directly in the external application

ATC³ Highest measuring accuracy
 'Absolute Temperature Calibration' for manual compensation of a temperature difference, 3-point calibration

Refrigeration Technology

Consistent cooling capacity
 Easily removable venting grid for quick and easy cleaning

ACC 100 % Cooling capacity
 'Active Cooling Control' for cooling available throughout the entire working temperature range, fast cool-down even at higher temperatures

Energy saving cooling
 Proportional cooling control for automatic adjustment of cooling power or temporary switch-off of compressor as needed to save up to 90 % energy in comparison to unregulated cooling machines

Technical Features

INTELLI PUMP Intelligent pump system
 Reliable and consistent pump capacity, electronically adjustable pump stages or pressure value, automatic adjustment of pump capacity to viscosity

ETHERNET Communication via networks
 For the remote control of instruments via Ethernet networks, full access to all functions of the unit via a networkcapable PC

USB Intelligent communication
 USB connection for data exchange (e.g. service data) or for wireless remote control via WirelessTEMP®

SmartCard Data exchange via SD-Card
 For data exchange (e.g. service data) via SD memory card

RS232 RS485 Connections according to standard
 RS232/RS485 dual-interface for serial data transmission according to EIA-485 industry standard (2-wire bus technology), upgradable with Profibus DP

8x60 Comfortable program control
 Integrated programmer for the execution of time and temperature dependant profiles, 8 temperature profiles with 60 steps max., with real time clock

Quiet as a whisper
 Efficient components produce only a minimal sound decibel level

Space-saving footprint
 All connections as well supply and exhaust air are located at the front or rear, no venting grids on the sides, units can be placed close to each other or the application

40° Continuous operation up to +40 °C
 Robust temperature control instrument, continuous operation even at ambient temperatures of up to +40 °C

EASY Transport Easy transport by one person
 Ergonomic design facilitates moving and positioning by one person

Filling level at a glance
 Backlit indicator for selected pump stages and filling volume

Warning & Safety Functions

Early warning system for high/low temperature limits
 Maximum safety for applications, optical and audible signal when limits are exceeded.

Duplicate safety
 Adjustable high temperature cut-off for internal tank and for integrated expansion vessel

S3 For flammable bath fluid
 Classification III (FL) according to DIN 12876-1

BLACK BOX Quick support
 If an error occurs, the integrated Black-Box function permits fast diagnosis by the JULABO service team