
OPERATING MANUAL

Ultra-Low Refrigerated Circulator

FP51-SL

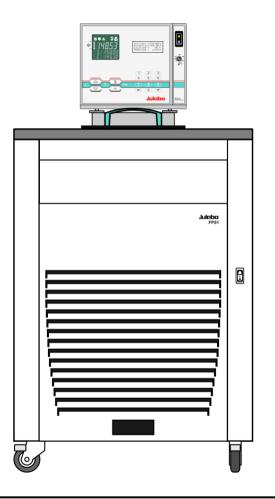



JULABO USA, Inc. 884 Marcon Boulevard Allentown, PA 18109 Phone: +1(610) 231-0250 Fax: +1(610) 231-0260 info@julabo.us www.julabo.us

1.950.2823.us.V08

01/22

Congratulations!

You have made an excellent choice.

JULABO thanks you for the trust you have placed in us.

This operating manual has been designed to help you gain an understanding of the operation and possible applications of our circulators. For optimal utilization of all functions, we recommend that you thoroughly study this manual prior to beginning operation.

The JULABO Quality Management System



Temperature control devices for research and industry are developed, produced, and distributed according to the requirements of ISO 9001 and ISO 14001. Certificate Registration No. 01 100044846

Unpacking and inspecting

Unpack the circulator and accessories and inspect them for possible transport damage. Damage should be reported to the responsible carrier, railway, or postal authority, and a damage report should be requested. These instructions must be followed fully for us to guarantee our full support of your claim for protecting against loss from concealed damage. The form required for filing such a claim will be provided by the carrier.

Important: keep operation manual for future use

TABLE OF CONTENTS

Operating manual	5
1. Intended use	5
1.1. Description	5
2. Operator responsibility – Safety recommendations	6
2.1. Disposal	7
2.2. Technical specifications	9
Operating instructions	11
3. Safety notes for the user	11
3.1. Explanation of safety notes	11
3.2. Explanation of other notes	11
3.3. Safety recommendations	12
4. Operating controls and functional elements	14
5. Preparations	18
5.1. Installation	18
5.2. Bath fluids	18
5.3. Temperature application to external systems	20
5.3.1.Tubing	21
5.4. Filling / draining	22
6. Operating procedures	23
6.1. Power connection	23
6.2. Switching on / Selecting the language	23
6.2.1.Switching on the circulator	23
6.2.2.Switching on the cooling machine	24
6.3. Start - Stop	24
7. Setting of temperatures	25
8. Safety installations, warning functions	25
8.0.1.Early warning system, low level protection	
9. Menu functions	
9.1. Start of a program	
9.1.1.Status at the end of a program	
9.1.2. Indication after successful start	
9.1.3.Editing after Start	
9.1.4.Interrupting a program	
9.1.5. Interruption after a power failure	
9.1.6.Termination of a program	
9.2. Program administration, creation	
9.3. Setting the pump pressure	
9.4. Configuration	
9.4.1.Remote control via the serial interface	
9.4.2.Keypad control or setpoint setting via the analog input	
9.4.3. Autostart	
9.4.4.Off-Mode	

9.4.5.ActVar - actuating variable	47
9.4.6.Setting of clock and date	48
9.4.7.Language	48
9.4.8.Unit	48
9.4.9.Program type	48
9.4.10. Reset	48
9.5. Control	49
9.5.1. Control internal / external	51
9.5.2.Selftuning	52
9.5.3.Dynamik intern	52
9.5.4.CoSpeed - extern	53
9.5.5.Control parameters – internal/external	53
9.6. Serial Interface	55
9.7. ATC Absolute Temperature Calibration, 3-point calibration	56
9.7.1.ATC Fühler - intern / extern	59
9.7.2.ATC Status - yes / no	59
9.7.3.Type 1 -/ 2 -/ 3 point	59
9.7.4.Example: 3-point calibration for internal control	59
9.8. Limits	62
9.8.1.Limits for internal control and external control	63
9.8.2.Limits for external control	64
9.9. Analog inputs/outputs	65
9.9.1.Control connector	69
9.9.2.Outputs of the connector - REG+E-PROG	70
9.9.3.Input of the connector - REG+E-PROG	71
9.9.4.External Stand-by input	73
9.9.5.Alarm-output	74
10. Troubleshooting guide / Error messages	75
11. Electrical connections	79
12. Remote control	82
12.1. Setup for remote control	82
12.2. Communication with a PC or a superordinated data system	82
12.3. List of commands	83
12.4. Status messages	87
12.5. Error messages	87
13. Installation of electronic module with analog connectors	
14. JULABO Service – Online remote diagnosis	89
15. Cleaning / repairing the unit	90
16. WARRANTY PROVISIONS	91

Operating manual

1. Intended use

JULABO circulators have been designed to control the temperature of specific fluids in a bath tank. The units feature pump connections for temperature control of external systems (loop circuit).



JULABO circulators are not suitable for direct temperature control of foods, semiluxury foods and tobacco, or pharmaceutical and medical products. Direct temperature control means unprotected contact of the object with the bath medium (bath fluid).

microprocessor technology allows to set and to store different values that can be indicated on the VFD COMFORT-DISPLAY and LCD DIALOG-DISPLAY. Three menu keys facilitate adjusting setpoints, warning and safety functions and menu

✓ The circulators are operated via the splash-proof keypad. The implemented

The integrated programmer allows storing and running temperature and time-

1.1. Description





The control electronics including "ICC - Intelligent Cascade Control" automatically adapts the heat supplied to the thermal requirements of the bath.
 The TCF - Temperature Control Features allow the user to have access to all

functions.

dependent processes.





mode and the chance to manually adjust or adapt control to the specific application.
 ✓ Absolute Temperature Calibration (ATC3) provides high temperature stability in the bath. With the 3-point calibration an offset is adjusted at three temperatures to ensure an accurate temperature pattern at the selected spot in the bath over the full temperature range.

important temperature control parameters. This means: Full control on the control



Pt100

SMART

PUMP

- Electrical connections:
 - The serial interface, switchable from RS232 to RS485, allows modern process technology without additional interface.
 - Connection for Pt100 external sensor for external temperature measurement and control.
 - Alarm output for external alarm message or control of JULABO refrigerating baths or solenoid valve (cooling water).
 - The electronic module (option) provides 3 further analog connections (alarm input, standby input, recorder output, programmer input).
- ✓ The excess temperature protection conforming to IEC 61010-2-010 is a safety installation independent from the control circuit. This protection can be indicated and set on the LCD DIALOG-DISPLAY.
- ☑ The early warning system for low level signals that bath fluid needs to be refilled before the low level protection conforming to IEC 61010-2-010 causes a complete shut-down of the main functional elements.
- ☑ Intelligent pump system: The pump capacity (electronically adjustable via the motor speed) enables to adapt to varying conditions for internal and external temperature applications.

2. Operator responsibility – Safety recommendations

The products of JULABO ensure safe operation when installed, operated, and maintained according to common safety regulations. This section explains the potential dangers that may arise when operating the circulator and also specifies the most important safety precautions to preclude these dangers as far as possible.

- > The operator is responsible for the qualification of the personnel operating the units.
- The personnel operating the units should be regularly instructed about the dangers involved with their job activities as well as measures to avert these dangers.
- Make sure all persons tasked with operating, installing, and maintaining the unit have read and understand the safety information and operating instructions.
- When using hazardous materials or materials that could become hazardous, the circulator may be operated only by persons who are absolutely familiar with these materials and the circulator. These persons must be fully aware of possible risks.

If you have any questions concerning the operation of your unit or the information in this manual, please contact us!

Contact	JULABO USA, Inc.	Phone:	+1(610) 231-0250
	884 Marcon Boulevard	Fax:	+1(610) 231-0260
	Allentown, PA 18109		info@julabo.us
			www.julabo.us

Safety instructions for the operator:

- You have received a product designed for industrial use. Nevertheless, avoid strikes to the housing, vibrations, damage to the operating-element panel (keypad, display), and contamination.
- Make sure the product is checked for proper condition regularly (depending on the conditions of use). Regularly check (at least every 2 years) the proper condition of the mandatory, warning, prohibition and safety labels.
- Make sure that the mains power supply has low impedance to avoid any negative effects on instruments being operated on the same mains.
- This unit is designed for operation in a controlled electromagnetic environment. This means that transmitting devices (e.g., cellular phones) should not be used in the immediate vicinity. Magnetic radiation may affect other devices with components sensitive to magnetic fields (e.g., monitors). We recommend maintaining a minimum distance of 1 m.
- > Permissible ambient temperature: max. 40 °C, min. 5 °C.
- > Permissible relative humidity: 50% (40 °C).
- > Do not store the unit in an aggressive atmosphere.
- > Protect the unit from contamination.
- > Do not expose the unit to sunlight.

Appropriate operation

Only qualified personnel is authorized to perform configuration, installation, maintenance and repairs of the circulator.

Routine operation can also be carried out by untrained personnel who should however be instructed by trained personnel.

Use:

The bath can be filled with flammable materials. Fire hazard!

There might be chemical dangers depending on the bath medium used.

Observe all warnings for the used materials (bath fluids) and the respective instructions (safety data sheets).


Insufficient ventilation may result in the formation of explosive mixtures. Only use the unit in well ventilated areas.

Only use recommended materials (bath fluids). Only use non-acid and non corroding materials.


When using hazardous materials or materials that could become hazardous, **the operator must** affix the enclosed safety labels to the front of the unit so they are highly visible:

If this unit is intended for use within the United States of America, all 2 warning labels **must** be affixed to the housing of the unit prior to use.

Directions for the positioning of the individual warning labels are enclosed with the warning labels included in the delivery. Warning labels must be easily visible to users.

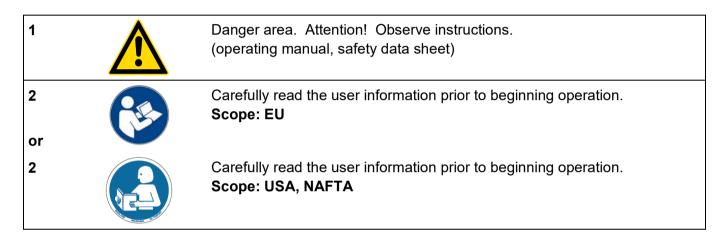

Particular care and attention is necessary because of the wide operating range. There are thermal dangers: Burn, scald, hot steam, hot parts and surfaces that can be touched.

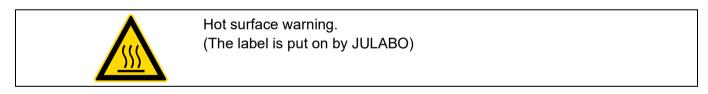

Observe the instructions in the manuals for instruments of a different make that you connect to the circulator, particularly the respective safety recommendations. Also observe the pin assignment of plugs and technical specifications of the products.

2.1. Disposal

The circulator contains a back-up battery that supplies voltage to memory chips when the unit is switched off. Do not dispose of the battery with household waste!

Depending on battery regulations in your country, you might be obliged to give back used or defect batteries to gathering places.

The product may be used with oil as bath fluid. These oils fully or partially consist of mineral oil or synthetic oil. For disposal, observe the instructions in the safety data sheets.

This unit contains refrigerants, which at this time are not considered harmful to the ozone layer. However, over the long operating period of the unit, disposal rules may change. Therefore, only qualified personnel should handle the disposal.



Contact an authorized waste management company in your country. Disposal with household waste (unsorted waste) or similar collections of municipal waste is not permitted!

2.2. Technical specifications

Performance specifications measured in accordance with DIN12876. Cooling capacities up to 20°C measured with ethanol; over 20°C with thermal oil unless specified otherwise. Performance specifications apply at an ambient temperature of 20°C. Performance values may differ with other bath fluids.

			-
Ultra-Low Refrigerated Circula	tor		FP51-SL
Working temperature range		°C	-51 200
Temperature stability		°C	±0,05
Cooling capacity		°C	<u>+20 0 -20 -40</u>
		kW	2.0 1.5 1.0 0.26
Cooling compressor			1-stage
Refrigerant			R507
Overall dimensions	(WxDxH)	cm	46x55x89
Bath opening	(WxL)	cm	18x12
Bath depth		cm	20
Filling volume		liters	7,5 12
Weight		kg	90
Mains power connection	z	V/ Hz	400 ±10 % / 50 / 3PNPE
Current draw	(at 400 V)	А	16
Mains power connection	(at 208 V)	V/ Hz	208 -5%; +10% 60 / 3PPE
Current draw	(at 230 V)	А	16
Mains power connection		V/ Hz	230 ±10 % / 50 / 3PPE
Current draw	(at 230 V)	А	16

			SL
Temperature selection			digital
via keypad			indication on LCD DIALOG-DISPLAY (°C/°F)
remote control via PC			indication on monitor
Temperature indication			VFD COMFORT-DISPLAY (°C/°F)
Resolution		°C	0.01
ATC3	INT / EXT	°C	±3 / ±9
Temperature control			ICC - Intelligent Cascade Control
Heater wattage (at 230 V	√)	kW	3.0
Electronically adj. pump	capacity	stages	1 4
Flow rate	at 0 bar	l/min	22 26
Max. pressure	at 0 liters	bar	0.7
Max. suction	at 0 liters	bar	0.4
Mains power connection		V/ Hz	208 ±10 % / 230 ±10 % / 50 / 60
Current draw	(at 208 V / 230 V)	А	13 / 14

Electrical connections:

External alarm device 24-0 V DC / max. 25 mA Computer interface RS232 or RS485 External Pt100 sensor

Optional for HL, SL

(Order No. 8900100 Electronic module with analog connections)			
Programmer input -1	00 °C to 400 °C = 0 -	10 V or 0 - 20 mA or 4 - 20 mA	
Input for the signal o	f a flow meter or exte	rnal manipulated variable	
Temperature recorder outputs $0 - 10 \vee (0 \vee = -100 \degree C, 10 \vee = 400 \degree C$			
	0 - 20 mA (0 mA = -	100 °C, 20 mA = 400 °C)	
	4 - 20 mA (4 mA = -	100 °C, 20 mA = 400 °C)	
Standby input	for external emerger	ncy switch-off	
Alarm output	for external alarm si	gnal	

Safety installations according to IEC 61010-2-010:

Excess temperature protection	adjustable from 0 °C 320 °C
Low liquid level protection	float switch
Classification according to DIN 12876-1	class III

Supplementary safety installations

Early warning system for low level	float switch
High temperature warning function	optical + audible (in intervals)
Low temperature warning function	optical + audible (in intervals)
Supervision of working sensor	plausibility control
Reciprocal sensor monitoring between	
working and safety sensors	difference >35 K
Alarm message	optical + audible (permanent)
Warning message	optical + audible (in intervals)

Environmental conditions according to IEC 61 010-1:

Use indoors only. Altitude up to 2000 m - normal zero. Ambient temperature: see Technical specifications Humidity: Max. relative humidity 80% for temperatures up to +31 °C, linear decrease down to 50% relative humidity at a temperature of +40 °C, max. permissible mains fluctuations, see Technical specifications.

Protection class according to IEC 60 529	IP21
The unit corresponds to Class	I
Overvoltage category	П
Pollution degree	2



Caution:

The unit is not suitable for use in explosive environment

EMC requirements

The device is an ISM device of group 1 per CISPR 11 (uses HF for internal purposes) and is classified in class A (industrial and commercial sector).

Notice:

- Devices of class A are intended for the use in an industrial electromagnetic environment.
- When operating in other electromagnetic environments, their electromagnetic compatibility may be impacted.
- This device is not intended for the use in living areas and cannot guarantee adequate protection of the radio reception in such environments.

Operating instructions

3. Safety notes for the user

3.1. Explanation of safety notes



In addition to the safety warnings listed, warnings are posted throughout the operating manual. These warnings are designated by an exclamation mark inside an equilateral triangle. "Warning of a dangerous situation (Attention! Please follow the documentation)."

The danger is classified using a signal word.

Read and follow these important instructions for averting dangers.



Warning:

Describes a **possibly** highly dangerous situation. If these instructions are not followed, serious injury and danger to life could result.



Caution:

Describes a **possibly** dangerous situation. If this is not avoided, slight or minor injuries could result. A warning of possible property damage may also be contained in the text.



Notice:

Describes a **possibly** harmful situation. If this is not avoided, the product or anything in its surroundings can be damaged.

3.2. Explanation of other notes

	Note! Draws attention to something special.
í	Important! Indicates usage tips and other useful information.
	This icon is used in the operating instructions to indicate flashing values or parameters which have to be set or confirmed.

3.3. Safety recommendations

Follow the safety instructions to avoid personal injury and property damage. Also, the valid safety instructions for workplaces must be followed.



- Only connect the unit to a power socket with an earthing contact (PE protective earth)!
- The power supply plug serves as a safe disconnecting device from the line and must always be easily accessible.
- Place the unit on an even surface on a base made of nonflammable material.
- Do not stay in the area below the unit.
- Make sure you read and understand all instructions and safety precautions listed in this manual before installing or operating your unit.
- Set the excess temperature safety installation at least 25 °C below the fire point of the bath fluid.
- Observe the limited working temperature range when using plastic bath tanks.
- Never operate the unit without bath fluid in the bath.
- Pay attention to the thermal expansion of bath oil during heating to avoid overflowing of the fluid.
- Prevent water from entering the hot bath oil.
- Do not drain the bath fluid while it is hot! Check the temperature of the bath fluid prior to draining (e.g., by switching the unit on for a short moment).
- Use suitable connecting tubing.
- Avoid sharp bends in the tubing, and maintain a sufficient distance from surrounding walls.
- Make sure that the tubing is securely attached.
- Regularly check the tubing for material defects (e.g., for cracks).
- Never operate damaged or leaking units.
- Always turn off the unit and disconnect the mains cable from the power source before performing any service or maintenance procedures, or before moving the unit.
- Always turn off the unit and disconnect the mains cable from the power source before cleaning the unit.
- Always empty the bath before moving the unit.
- Transport the unit with care.
- Sudden jolts or drops may cause damage in the interior of the unit.
- Observe all warning labels.
- Never remove warning labels.
- Never operate units with damaged mains power cables.
- Repairs are to be carried out only by qualified service personnel.



Some parts of the bath tank and the pump connections may become extremely hot during continuous operation. Therefore, exercise particular caution when touching these parts.



Caution:

The temperature controlling i.e. of fluids in a reactor constitutes normal circulator practice.

We do not know which substances are contained within these vessels. Many substances are:

- inflammable, easily ignited or explosive
- hazardous to health
- environmentally unsafe

i.e.: dangerous

The user alone is responsible for the handling of these substances! The following questions shall help to recognize possible dangers and to reduce the risks to a minimum.

- Are all tubes and electrical cables connected and installed? Note:
 - sharp edges, hot surfaces in operation, moving machine parts, etc.
- Do dangerous steams or gases arise when heating? Is an exhaust needed when working?
- What to do when a dangerous substance was spilled on or in the unit? Before starting to work, obtain information concerning the substance and determine the method of decontamination.



Notice:

Check the safety installations at least twice a year!

- Excess temperature protection according to IEC 61010-2-010. With a screwdriver turn back the adjustable excess temperature protection until the shut-down point (actual temperature).
- Low level protection according to IEC 61010-2-010. To check the function of the float, it can be manually lowered with a screwdriver for example.





WARNING


This product contains chemicals known to the state of California to cause cancer, birth defects or other reproductive harm.

4. Operating controls and functional elements

Front view

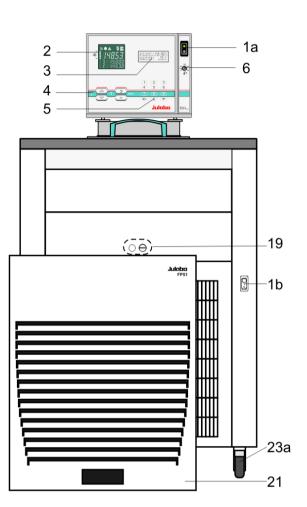

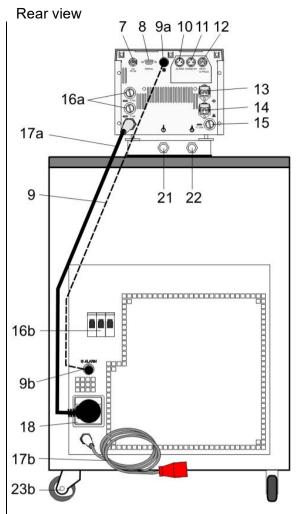

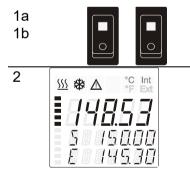

Mains power switch, illuminatedfor circulatorMains power switch, illuminatedfor cooling machine

VFD COMFORT-DISPLAY

Header: Control indicators

- Line 1: Actual value internal or external The display is depending on the selected control mode in the menu > Control < (internal or external).
- Line 2: Working temp. setpoint, constantly S xxx.xx
- Line 3: Actual value (E = external or I = internal) Alternating with the display in line 1

Use the keys 🔘 to indicate further values in line 3


PI Capacity in % - with manipulated variable set to >control<*

PS Capacity in % - with manipulated variable set to >serial<* or >eprog<*

- H Heater capacity in Watts
- U Mains voltage Volts
- F Flow rate in liters/minute

(providing EPROG input set to >Flow rate<)

*refer to >MENU / CONFIGIGURATIO> page 32

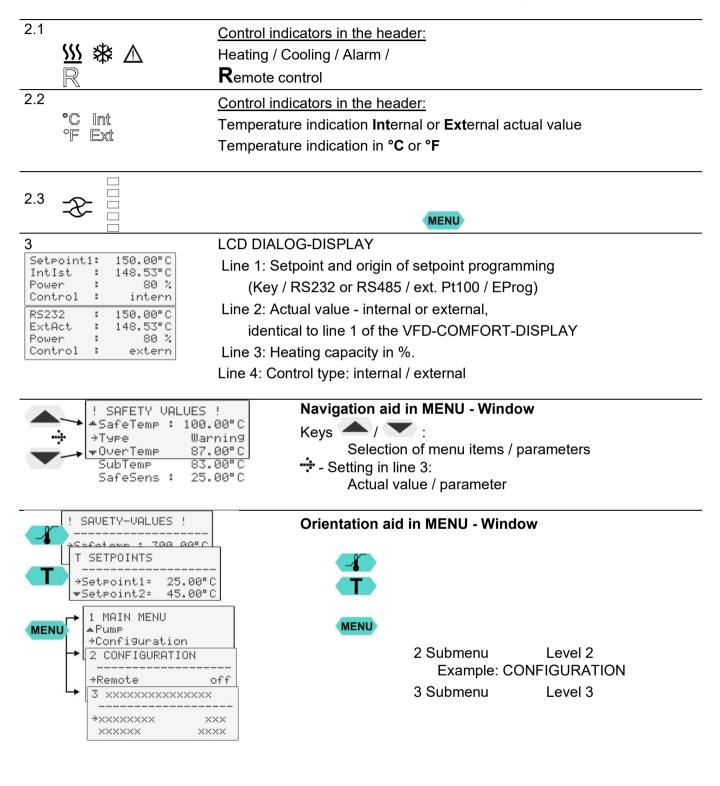

4	Keypad 1	Navigation keys
1.1	ок	1. Key: >OK< Start / Stop (pump / heater)
		2. >OK< in the menu Menu item / select submenu for setting
		Save set value
		Save selected parameter
	~	A beep signals the end of setting
		After the actions Start, Stop and change from VFD Display to standard
		display the key OK is locked for a short time.
4.2	•	The above graph "front side" shows an example for standard display.1. Key: >ReturnStop (pump / heater)
4.2	J	Special display on VFD Display leave line 3.
		2. >Return< in the menu one menu level down
		Correction function for parameters or values (prior to OK)
		• (D)
		immediately back to standard display
	r and the second s	οκ (P) - () (P) icon for "keep key pressed down".
		the pressed down.
4.3		1. Key: >Up / Down <temperature decrease="" increase="" setpoint<="" td="" –=""></temperature>
-		Push key quickly for single steps,
		Keep key pressed for fast change.
		2. >Up/Down< in the menu selection of menu items / parameters
		Menu keys
4.4		Key: start the menu > warning and safety values<
4.5	T	Key: start the menu >temperature setpoints<
4.6	MENU	Key: display of MENU structure
5	Keypad 2	
•		Numerie keyned: numerele 0 te 0
	0.9	Numeric keypad: numerals 0 to 9
	-	minus / decimal point
		Backspace key SetPoint1: 155 °C
		on LCD DIALOG-DISPLAY.
6	⁸⁰ ⁴⁰ ⁴⁰ ⁵ ³²⁰ ⁵ ⁵	Adjustable excess temperature protection according to IEC 61010-2-010

Rear view

7	ext Pt100	Socket for external measurement and control sensor or external setpoint programming
8	ہ ‱ ہ SERIAL	Interface RS232 / RS485: remote control via personal computer
9		Control cable of JULABO refrigerated circulator
9a	*	Socket: output for alarm messages
9b	•	Socket:

Optio	n: Electronic modu	le Order No. 8 900 100		
		The circulator automatically recognizes the connected electronic module.		
10		Alarm output (for external alarm signal)		
	ALARM			
11		Standby input (for external emergency switch-off)		
10	STAND-BY			
12		Programmer input and temperature recorder output		
	REG+E-PROG			
13	€	Connector for supplementary pump or solenoid valve 230 V / max. 1.25 A		
14		No control voltage in the -OFF- condition Connector solenoid valve		
14		1. No control voltage in the -OFF- condition		
15		2. Configurable (Refer to page 65)		
15		Fuses (for connectors 13, 14), T1.25A		
16		Mains fuses, T16A		
16b		Mains circuit breakers (resettable) 16 A		
17a		Mains power cable with plug for circulator		
17b		Mains power cable with plug cooling machine		
18	\odot	Built-in mains outlet for connection of circulator (230 V / 16 A)		
19	$\bigcirc \Theta \bigcirc \blacksquare$	Drain screw with drain connection		
20		Venting grid, removable		
21 22		Pump connections U suction pump O pressure pump		
23a		Castor with brake (at the front)		
23b		Castor without brake (at the back)		

5. Preparations

5.1. Installation

- Place the unit on an even surface on a pad made of **non-flammable** material. Using the castors move the unit to the intended location. For better stability, apply the holding breaks on the front casters.
- Cooling machine, pump motor and electronics produce intrinsic heat that is dissipated via the venting openings! Never cover these openings!
- Keep at least 20 cm of open space on the front and rear venting grids.
- The place of installation should be large enough and provide sufficient air ventilation to ensure the room does not warm up excessively because of the heat the instrument rejects to the environment. (Max. permissible ambient temperature: 40 °C).
 For a fault (leakage) in the refrigeration system, the standard EN 378 prescribes a certain room space to be available for each kg of refrigerant.
 The refrigerant quantity is specified on the type plate.
 > For 0.49 kg of refrigerant R507, 1 m³ of space is required.
- Do not set up the unit in the immediate vicinity of heat sources and do not expose to sun light
- Before operating the unit after transport, <u>wait about one hour after</u> <u>setting it up</u>. This will allow any oil that has accumulated laterally during transport to flow back down thus ensuring maximum cooling performance of the compressor.

5.2. Bath fluids



Caution:

Carefully read the safety data sheet of the bath fluid used, particularly with regard to the fire point!

If a bath fluid with a fire point of \leq 65 °C is used, only supervised operation is possible.

Water: The quality of water depends on local conditions.

- Due to the high concentration of lime, hard water is not suitable for temperature control because it leads to calcification in the bath.
- Ferrous water can cause corrosion even on stainless steel.
- Chloric water can cause pitting corrosion.
- Distilled and deionized water is unsuitable. Their special properties cause corrosion in the bath, even in stainless steel.

Recommended bath fluids:

Bath fluid	Temperature range
soft/decalcified water	5 °C to 80 °C
mixture water/glycol, mixture 1:1	-20°C to 50°C

JULABO bath fluids

JULABO		Thermal	Thermal	Thermal
Description		G	HY	H5
Order Number	10 liters	8 940 124	8 940 104	8 940 106
	5 liters	8 940 125	8 940 105	8 940 107
Temperature range	°C	-30 80	-80 55	-50105
Flash point	°C		78	124
Fire point	°C		80	142
Color		light yellow	clear	clear

JULABO		Thermal	Thermal
Description		H10	H20S
Order Number	10 liters	8 940 114	8 940 108
	5 liters	8 940 115	8 940 109
Temperature range	°C	-20 180	0 220
Flash point	°C	190	230
Fire point	°C	216	274
Color		clear	light brown



See website for list of recommended bath fluids.

ATTENTION:

The maximum permissible viscosity is 70 mm² /s·



Caution:

Fire or other dangers when using bath fluids that are not recommended:

Use only non-acidic and non-corroding bath fluids. JULABO assumes no liability for damage caused by the selection of an unsuitable bath liquid.

Unsuitable bath fluids are fluids which, e.g.,

- are highly viscous (much higher than recommended at the respective working temperature)
- have a low viscosity and have creep characteristics
- have corrosive characteristics or
- tend to crack.
- No liability for use of other bath fluids!


5.3. Temperature application to external systems



Caution: Securely attach all tubing to prevent slipping.

If the circulator is operated without external system, close the pump connector (22) with the cap nut.


Temperature application to external, closed systems

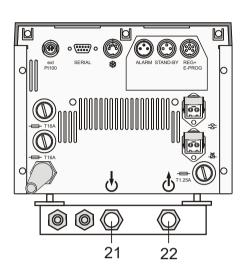

The circulator is used for temperature application to external, closed systems (loop circuit) with simultaneous temperature application in the circulator bath.

- Unscrew the M16x1 collar nuts on the pump connectors with a 19 mm (3/4") wrench and remove the sealing disks. Using the collar nuts, screw on the tubing connection fittings (for tubing 8 mm or 12 mm in diameter) delivered with the unit and tighten firmly. (Pressure pump: 22, suction pump: 21)
- Push on the tubing, and secure with tube clamps.
- Attach the tubing to the connectors of the external closed system, e.g., an instrument with a pressure-resistant temperature jacket or a temperature coil, and fasten with tube clamps to prevent slipping.

Temperature application to external, open systems

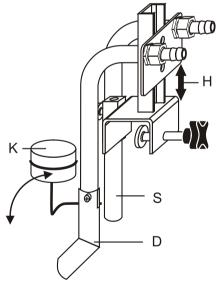

- S = Suction pump connection
- D = Pressure pump connection
- K = Float
- H = Height adjustment

The circulator is equipped with both a pressure and suction pump for external temperature application in open systems.


Differing flow rates of the pressure and suction pumps should be compensated. To maintain a constant liquid level, the JULABO "D+S" Level Adapter is recommended for the external bath tank. The flow rate of the pressure pump will be then regulated by a built-in float device. The liquid level may be changed by a height adjustment on the "D+S" Level Adapter.

Accessory: "D+S" Level Adapter Order No. 8 970 410

Important:

- The liquid level should be equal in the internal and external baths (absolute height).
- (i) If you take out samples (for example Erlenmeyer flasks) from the external bath, turn the circulator off with the Start/Stop key.

Backflow safety device

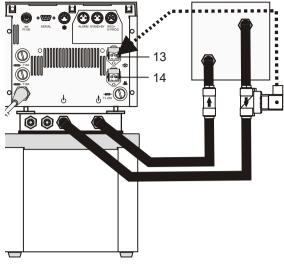

2 INPUTS/OUTPUTS			
→SV-OUT Backflow			
▼Chani ActInt			
(see page 65			

5.3.1. Tubing

Recommended tubing:

If the liquid levels in the circulator bath and the external system are at different heights, overflowing must be prevented after the power has been turned off.

Solution Flood hazard!

For this reason, solenoid valves for loop circuit or shut-off valves can be integrated in the loop circuit.

Connect the valve to the connector (13).

 If socket (14) is used: In menu >Inputs/Outputs< set the menu item >SV-Out< to >Backflow<.

Order No. Description

- 8 980 701 Set of solenoid valves for loop circuit up to +100 °C
- 8 970 456 Shut-off valve (suitable up to +90 °C)
- 8 970 457 Shut-off valve (suitable up to +200 °C)

ige
°C
С
•



Warning: Tubing:

At high working temperatures the tubing used for temperature application and cooling water supply represents a danger source.

A damaged tubing line may cause hot bath fluid to be pumped out within a short time. This may result in:

- Burning of skin
- Difficulties in breathing due to hot atmosphere

Safety recommendations

- Employ suitable connecting tubing.
- Make sure that the tubing is securely attached.


- Avoid sharp bends in the tubing, and maintain a sufficient distance from surrounding walls.
- Regularly check the tubing for material defects (e.g. for cracks).
- Preventive maintenance: Replace the tubing from time to time.

5.4. Filling / draining



Notice:

- Pay attention to the thermal expansion of bath oil during heating to avoid overflowing of the liquid.
- Do not drain the bath fluid while it is hot! Recommendation: Temperature range 5 °C to 40 °C Check the temperature of the bath fluid prior to draining (by switching the unit on for a short moment, for example).
- Store and dispose the used bath fluid according to the laws for environmental protection.

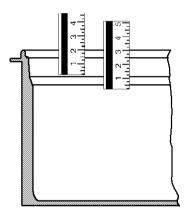


Filling

Take care that no liquid enters the interior of the circulator.

- Recommended maximum filling level with water as bath fluid: 30 mm below the tank rim
- Recommended maximum filling level with bath oils: 40 mm below the tank rim
- (1) After filling, immerse the samples in the bath or place the lid on the bath, in case the opening is not to be used.
- The circulator provides an early warning system for low level (description – please refer to page 29) that may be triggered when changing samples in the bath.



- Turn off the unit and disconnect the mains cable from the power source.
- Hold the venting grid, pull out and remove.
- Slide a short piece of tube onto the drain connection and hold it into a container.
- Loosen the drain screw a few turns and drain the unit completely.
- Tighten the drain tap.

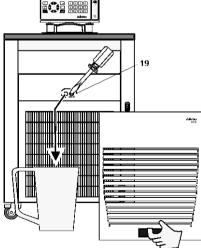

6. **Operating procedures**

6.1. **Power connection**


Caution:

- Only connect the unit to a power socket with earthing contact (PE protective earth)!
- The power supply plug serves as safe disconnecting device from the line and must • be always easily accessible.
- Never operate equipment with damaged mains power cables.
- Regularly check the mains power cables for material defects (e.g. for cracks).
- We disclaim all liability for damage caused by incorrect line voltages!



Check to make sure that the line voltage matches the supply voltage specified on the identification plate.

Mains power connection:

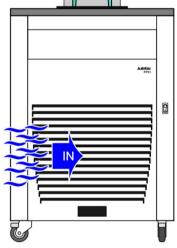

3x400V/N/PE

V/N/PE

400

- 3x230V/PE - 1 L1 = 230 V/PE L2 = 12 -L3 ⊏ 13 5 N PE _____gngb / gnye PE _____angb / anvel
- (i) Note the direction of rotation of the built in fan! For proper cooling of the compressor, the air must be drawn in from the front of the unit.
- Connect the circulator with mains power cable (17a) to the mains • outlet (18).
- Connect the control cable (9) to the sockets (9a, 9b).
- Connect the refrigerated circulator with mains power cable (17b) to the mains socket.

Switching on / Selecting the language 6.2.

6.2.1. Switching on the circulator



JULABO 111 U1.xx xxxx 150.00°C Setpoint1: 148.53°C IntIst 5 Power 2 80 % Control ÷ intern R 50.00

T

24.64

- Switching on:
- Turn on the mains power switch (1a).

(i) The unit performs a self-test.

Then the software version (example: V 1.xx-xxxx) appears. The display "OFF" or "R OFF" indicates the unit is ready to operate.

(i) The circulator enters the operating mode activated before switching the circulator off:

keypad control mode (manual operation) or remote control mode (operation via personal computer).

23

2 KONFIGURAT	ION
≜ Zeit/Datum →Sprache	deutsch
∀ Einheit	°C
2 CONFIGURAT <i>*</i> Time/Date	IUN

Selecting the language:

There are two options for the language of the LCD DIALOG-DISPLAY: German or English. Select the desired language in the menu >Configuration< under the submenu >Language/Sprache<.

Press the respective keys in the following order:



ок

6х up to submenu >Language/Sprache< ок

K deutsch / english Жок Э (У)

6.2.2. Switching on the cooling machine

Switching on:

• Switch on the cooling machine using the switch (1b).

(i) Control of the cooling machine:

With the mains switch (1b) turned on, the circulator automatically switches the cooling machine off and on.

- It is switched off if:
 - the current working temperature is increased by > 30 °C (cooling is not required).

- the heater operates at full power (> 800 W) for longer than 5 minutes.

It is switched on if:

- cooling is necessary to maintain the bath temperature. After switch-off, the cooling machine automatically switches on only after a delay of 5 minutes to protect the cooling compressor.

To save energy, turn off the cooling machine with the mains switch (1b) whenever cooling is not required.

6.3. Start - Stop



Start:

Press **OK** key. The actual bath temperature is displayed on the VFD COMFORT-DISPLAY. The circulating pump starts with a slight delay.

Stop:

Press OK key.

or

Keep () () key pressed.

The VFD COMFORT-DISPLAY indicates the message "OFF".



Press the **T** key to call up the menu for temperature selection.

3 different working temperatures can be adjusted. Their values are freely selectable within the operating temperature range.

① The temperatures can be set in start or stop mode.

Setting of working temperature in the

- 1. Press the key \frown . The value flashes \mathcal{W}
- 2. Select SETPOINT 1 or 2 or 3 using the key 💙 or 📥
- 3. Confirm by pressing the **OK** key.
- (1) The circulator uses the new working temperature value for temperature control.

Example: Adjustment/modification of the pre-setting of "SETPOINT 3"

- If the active Setpoint is changed, the new value is immediately used for the control of the working temperature. The heater control indicator flashes.
- Setpoint3: Change the value from 70.00°C to 80.00°C. See standard display line 1
- Use the numeric keypad to enter 8 and 5. Then press enter **OK** to store the value.

Notice: Refer to SETPOINT MAX / MIN in chapter "LIMITS"



Check the safety installations at least twice a year! Refer to page 13.

Settings for the excess temperature protection > SafeTemp< and for the warning functions > OverTemp< and > SubTemp< are made in a menu which is called up by pressing the key

Menu item > Type< allows choosing between a warning and an alarm cut-off for the menu items > OverTemp< and > SubTemp<.

Factory settings:

i a	iciory setting	<u>j</u> 3.			
Т	SETPOINTS				
·					
l⇒ś	Setpoint1	=	25.	00°	С
w S	Setpoint2	=	37.	00°	С
	Setpoint3	=	70.	00°	С

Setpoint3:	70.00°C
ActIst :	24.64°C
Power :	80 %
Control :	internal
Setpoint3:	85.∎0°C
ActIst :	24.64°C
Power :	80 %
Control :	internal




Warning:

The excess temperature protection **must** be set at least 25 °C below the fire point of the bath fluid used! In case of wrong setting there is a fire hazard!


We disclaim all liability for damage caused by wrong settings!

J

! SAFETY VALUES ! →SafeTemp : 100.00° wType Warni OverTemp 87.00° SubTemp 83.00° SafeSens : 25.00°	n⊈ ∕ Alarm temperature scale. C			
	 Press the key and by pressing the key or select 			
	the menu item \Rightarrow and confirm the by pressing the key \circ .			
	For setting proceed as follows.			
SafeTemp:	 Set the new cut-out value within 30 seconds with a screw driver. Exact display and setting on LCD display Recommendation: Set over temperature-protective setting >SafeTemp< 5 °C to 10 °C above the set point of the working temperature. 			
Туре:	Switch-over from warning to shutdown function			
	 Change the parameter by pressing the key and confirm by pressing the key OK . or pressing the key if the parameter is to be retained. Possible parameters: warning or alarm 			
OverTemp :	• Set value on numeric keypad and confirm by pressing OK			
SubTemp:	Example: 83.00 °C			
	Press the keys 8 and 3 and confirm by pressing OK			
	Indicated is the temperature value of the safety sensor.			
	SafeTemperature:			
	Setting range: 20 °C 320 °C			
	Indicated is the cut-out value set with a screwdriver on the excess temperature protection device.			
	This safety installation is independent of the control circuit. When the temperature of the bath fluid has reached the safety temperature, a complete shutdown of the heater and pump is effected. The alarm is indicated by optical and audible signals (continuous tone). The following error message appears on the VFD COMFORT-DISPLAY:			
26				

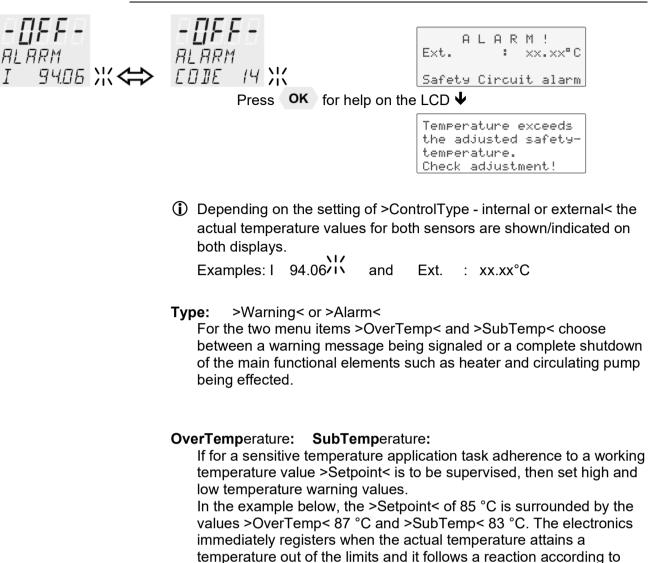

what is set in the menu item >-Type<.

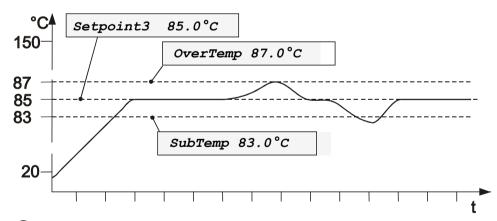

The warning functions are only triggered when the actual bath temperature, after start from the "OFF" or "rOFF" mode, lies within the set limits for 3 seconds.



A mere warning function with optical and audible warning signal (interval tone) A message appears on the VFD COMFORT-DISPLAY:



	1
VARNINE CODE DE	
OverTemp	

NABNING EDIE DY SubTemp

The LCD – DISPLAY shows the message:

Setpoint1: 150.00°C	Setpoint1: 150.00°C
IntAct : 148.53°C	
excess temperature!	low temperature!
-check limits-	-check limits -

Setting >ALARM< •

Temperature limit with shutdown of heater and circulating pump. An audible alarm sounds (continuous tone) and a message appears on the VFD COMFORT-DISPLAY:

- [] F F -	
ALARM COIE OB	
OverTemp	or

- [] F F -RLARM EDIE DY SubTemp

The LCD -DISPLAY shows the message

ALARM! Ext. : xx.xx*C	ALARM! Ext. : xx.xx*C	
Excess Temperature	Low Temperature!	
Press OK for help on the LCD		
	Ostrin 1. A sum su struct	

Actual temperature obove high temperature value. Check limits!

Actual temperature below low temperature value. Check limits!



8.0.1. Early warning system, low level protection




DBGM: 20306059.8


This low level protection is independent of the control circuit and is divided into two sections:

1. Switch in stage 1 recognizes a critical fluid level $\textcircled{\odot}$.

An audible warning (interval tone) sounds and a message appears on the VFD COMFORT-DISPLAY.




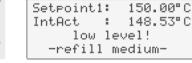

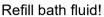

- Switch in stage 2 recognizes a low fluid level .
 If stage 2 of the low level protection device (according to IEC 61010-2-010) is triggered, a complete shutdown of the heater and circulating pump is effected.

A continuous alarm tone sounds and a message >ALARM< >CODE 01< appears on the VFD COMFORT-DISPLAY.

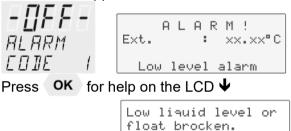

medium!

Turn off the unit with the mains switch, refill bath fluid and turn the unit on again!

Check float and fill

- 3. Float
- 4. Circulating pump
- 5. Heater



Warning:

For refill always use the same bath fluid type that is already in the bath. Bath oils must not contain any water contaminants and should be pre-heated to the actual bath temperature! Explosion hazard at higher temperatures!

MENU





1 MAIN MENU →Start Pro9ram ▼Mana9e Pro9rams Pump **Configuration** Control Interface serial ATCalibration Limits Inputs/Outputs

Menu level 1

1 MAIN MENU →Start Pro9ram ▼Mana9e Pro9rams	Start program This menu will start a previously set program.	Page 32
1 MAIN MENU ▲Start Program →Manage Programs ▼Pump	Administration and creation of programs	Page 37
1 MAIN MENU ▲Manage Programs →Pump ▼Configuration	Electronically adjustable pump capacity The capacity of the circulating pump is set by adjusting the m	Page 41 otor speed.
1 MAIN MENU ▲Pump →Configuration ▼Control	 Configuration of the unit Remote – on / off (remote control via RS232) Setpoint ext - Switch over external setpoint setting Autostart – AUTOSTART on / off Off-Mode – Pump on / off ACTVAR - Switch over input variable Time / Date – setting time and date Language – english / deutsch Unit – Temperature values can be displayed in °C or °F Program Type – Time or Gradient (grad) Reset – factory settings 	Page 42
1 MAIN MENU ▲Configuration →Control ↓Interface serial	Control characteristics and parameters ControlType - Control internal or external Selftuning Dynamic - internal CoSpeed – external Control parameters - XP-, TN-, TV- internal Control parameters - XP-, TN-, TV- XPU-, external	Page 49

The term "menu functions" refers to adjustments such as

1 MAIN MENU ▲Control →Interface serial ₩ATCalibration	Adjustable interface parameters BAUDRATE, PARITY, HANDSHAKE	Page 55
1 MAIN MENU ▲Interface serial →ATCalibration ↓Limits	ATC - Absolute Temperature Calibration, Sensor calibration internal sensor, Sensor calibration external sensor 3-point calibration	Page 56
1 MAIN MENU ATCalibration +Limits FINPUTS/OutPuts	Limitations of temperature and capacity Setpoint Max / Min - Maximum and minimum setpoint Heating Max – Set maximum heating Cooling Max – Set maximum cooling Internal Max / Min – Limitation of the temperature range Upper band / Lower band – Band limit	Page 63
1 MAIN MENU ▲Limits →Inputs/Outputs	Analog inputs/outputs Recorder output – CHANNEL 1, 2, 3 EPROG – External programmer input Ext-StBy - Stand-By input ALARM - output	Page 65

9.1. Start of a program

2 START PROGR	:AM
→Program Numb	er 0
▼at Step	0
Runs	1
Pro9ram End	StdBy
Run	now

The start menu of the integrated programmer allows one of six previously stored temperature programs to be called up and started. The profiles are started manually or via the integrated timer.

(i) Requirements:

- 1. Create a program. (refer to next chapter)
- 2. Return to the Start-MENU and confirm the desired setting of each MENU item with the key **OK**
- 3. Set a start time (>TIME< >DATE< >YEAR<) if the program is to be started by the internal timer.

>Program Number< Select the program that should be started. 0 5

	05	
> at Step<	Select the Step at which the program should be started. 0 59	
> Runs <	Select the number of program repetitions. 1 99	
> Program End<	Status at program end .	
	StdBy, PSetp, Setp 1, Setp 2, Setp 3	
> RUN <	Start time (now / time)	
Press the	key if a parameter is to be retained. (prior to OK)	

Level 2	Parameter /Value	
2 START PROGRAM →Program Number 0 ▼at Step 0	2 START PROGRAM Program Number 0	 Set a parameter using the numeric keypad and confirm by pressing OK. Possible parameters: 05
2 START PROGRAM AProgram Number 0 →at Step 0 ▼Runs 1 → or	2 START PROGRAM at Step 0	 Set a parameter using the numeric keypad and confirm by pressing OK Possible parameters: 0 59
2 START PROGRAM At Step 0 Aruns 1 Program End StdBy	2 START PROGRAM 	 Set a parameter using the numeric keypad and confirm by pressing OK Possible parameters: 1 99
2 START PROGRAM ARuns 1 →Program End StdBy TRUN now	2 START PROGRAM →Program End ♣StdBy	 Set desired parameters with and ок. Possible parameters: StdBy , PSetp, Setp 1, Setp 2 , Setp 3

1	M	IENU	$\mathbf{\Psi}$	
	1	MAIN	MENU	
			Program e Programs	

ок 🕁

	Parameter /Value	
2 START PROGRAM [♠] Program End StdB →RUN no 	IZ SIHKI PRUGRHM	 Set desired parameters with and OK. Confirm >now< with the OK key and the program will start immediately or start at the set time under parameter (time). Set time in the example below: 09. August 2009, 11:15 hrs
3 TIMER →hour:min 11:1 *Day.Mon 09.0 Year 200 RUN	в	Set the time for the start of the program in the submenu >3 TIMER<.
3 TIMER >hour:min 11:1 *Day.Mon 09.0		 Set starting time using the numeric keypad and confirm by pressing ОК. Example: (hh:mm) 1 1 5 + ОК
3 TIMER ≜hour:min 11:1 ⇒Day.Mon 09.0 ▼Year 200	B 3 1111EK	 Set day and month using the numeric keypad and confirm by pressing OK 0 9 - 0 8 OK
3 TIMER ★Day.Mon 09.0 →Year 200 ▼RUN	I S LIMER	 Set year using the numeric keypad and confirm by pressing OK. Example: (JJJJ) 2 0 0 9 + OK
3 TIMER ≜Year 200 →RUN	9 3 TIMER 	Confirm >Starten/Start?< by pressing ок and the program starts at the set time.
Starttime: 11:1 Date : 09.0	Display of waiting time unt The unit switches to waiting	il start: g mode and a flashing line "wait" appears on Y. The start time and current time are

***** Wait ***** Act. Time: 10:34:57

permanently indicated on the display.

It might be necessary to check the correct setting of the internal real time clock

(see MENU configuration)

9.1.1. Status at the end of a program



Setpoint	3:	80.00°C
IntAct	:	80.00°C
Power	:	3%
Control	:	internal

T SETPOINTS	5	
▲Setpoint1	=	25.00°C
⇒Setpoint2	=	58.00°C
▼Setpoint3	=	85.00°C
Setpoint2:		58.00°C
IntAct :		60.00°C
Power :		0%
Control :		intern



Program end StdBy , PSetp, Setp 1, Setp 2 , Setp 3

This parameter is used to determine if the unit switches to the -OFFcondition at the end of a profile or if temperature control is continued, and the working temperature setpoint that is to be used if temperature control is continued.

StdBy – If the end of the profile is reached, the circulator switches to the -OFF- condition. The VFD display indicates the working temperature setpoint that was displayed in line 2 before the program was started (example: S 25.00 °C).

PSetp –(Setpoint of the **P**ro**G**rammer) At the end of the program, the setpoint of the last program step is written to the position of setpoint 3 (example: 80.00 °C) and indicated in line 1 of the LCD display. The circulator operates and controls to this temperature until the key **OK** is pressed or a new profile is started.

Setp 1, Setp 2, Setp 3 – ("Setp" short for Setpoint)

Before starting the programmer, a temperature value is set for one of the three setpoints in the **menu**.

(example: setpoint 2 58.00 °C)

The circulator operates and controls to this temperature at the end of the profile until the key **OK** is pressed or a new program is started.

9.1.2. Indication after successful start

The started programmer displays the currently calculated setpoint in line 1. The value changes until the final temperature of the section is reached (temperature ramp).

1st line: Setpoint of the programmer

2 nd line:	Actual temperature value	
	for internal control = IntAct:	XXX.XX
	for external control = ExtAct:	XXX.XX
3 rd line:	Selected profile and the curren	t section (step

- 3^{rd} line: Selected profile and the current section (step). 4^{th} line: With the \checkmark \bigstar keys, the following information can be
- displayed in line 4.
 - Remaining time of the current section (standard display)
 - Total remaining time: profile x number of repetitions (All: Days/hours:minutes:seconds)

Particularities :

- If a total time of more than 999 days is calculated, only the hours are displayed by the unit.
 (All: hours:minutes:seconds)
- 2. If the time period of a section is edited <u>after</u> the start, then the remaining total time can no longer be displayed.

9.1.3. Editing after Start

The programmer allows a running program to be edited.

Examples:

Set	:	XX.XX°C
IntAct	:	XX.XX°C
ProgNr.	:2	Step: 2
Remain	:	03:32:25

Set	:	XX.XX°C
IntAct	8	XX.XX°C
ProgNr.	:	2 Step: 2
All	8	ttt/hh:mm:ss

3	Εľ	DITOR	>Pr9Nr.2
SP	1	(C xxx.xxC
ΤI	1		4 ××.××4
ST	:		# 0

Initiation and exit of the editing function:

After the start enter the submenu "3 EDITOR".by pressing the key MENU

See chapter "program administration, creation" page 37


Exit the editing modus anytime by pressing the key 🕥 .

Particularity for changes in a section that is currently in process.

- (i) If the target temperature is changed, the slope of the ramp is automatically calculated based on the remaining time of the section.
- (i) If the time period is changed, the slope of the ramp is automatically calculated based on the new total time of the section.
- (i) The "remaining total time" (profile x number of repetitions) as described on page 34 can no longer be displayed.

9.1.4. Interrupting a program

Set	:			Χ	Χ.	XX°	С
IntAct	8			Х	Χ.	XX°	C
ProgNr.	8	2	S	t	eP	:	2
***		Paus	e		**	*	

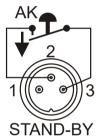

Set	:	XX.XX°	C
IntAct	:	XX.XX°	C
ProgNr	.:2	Step:	2
****	Star	ndBy ****	*

() Press the () key to interrupt or continue a profile.

The setpoint and time period set for the corresponding section are thus stopped at the values presently achieved.

The instrument is put on hold and the message "pause" flashes on the LCD DIALOG DISPLAY.

() A program can be interrupted or restarted by an external shut-off.

Important:

For this, in the >Inputs/Outputs < menu, the menu item >Ext. StBy < must be set to >active<

and in the >Configuration < menu, >Autostart < needs to be set

(i) Setpoint control and timer are interrupted by breaking the contact "AK".


The instrument is put on hold and the message "Stand-By" flashes on the LCD DIALOG-DISPLAY.

Notice: This is not an actual emergency switch-off.



Warning:

Following a power interruption, it would be possible in this condition for the instrument to restart automatically. The safety and warning functions of the instrument should always be used to their fullest capacity. See Warning page 46.

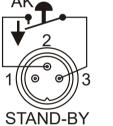

Starttime:

. ***** wait *****

Act. Time: 10:34:57

Date

9.1.5. Interruption after a power failure

11:15

09.08

It is possible to operate the programmer safely in case of a power failure.

Important:

For this, >Autostart< needs to be set in the >Configuration< menu.

- (i) If the AUTOSTART function is activated, the programmer starts again at a point approx. 20 seconds before the interruption took place. However, an uncontrolled change of the bath temperature has occurred.
- (i) Power failure when the instrument is on hold if start is effected via the built-in timer:
 - If the starting time is not yet reached, the programmer is put on hold again.

If the starting time is exceeded, the programmer starts immediately.

9.1.6. Termination of a program

Set	8	XX.XX°C
IntAct	8	XX.XX°C
ProgNr.	:2	Step: 2
Remain	8	03:32:25

2 START PROGRAM	
→Program Number	0
‴at Step	0
Runs	1
Program End 9	StdBy
Run	now

(1) A program can be terminated by pressing $\bigcirc \bigcirc$. The programmer switches back to the Start menu.

Further procedures are defined in the MENU >START PROGRAM<, menu item >Program End <.

Example: >Program End - StdBy<

The VFD COMFORT-DISPLAY displays "OFF".

Interruption after a power failure

In case of a power failure, the reaction of the circulator is the same as when switched off and on again with the mains switch.

The VFD COMFORT-DISPLAY displays "OFF".

Program administration, creation 9.2.

2	MANAGE	PROGRAMS
-		
⇒E	Edit	Pr9Nr.
T)elete	Pr9Nr.
F	^o rint	Pr9Nr.

The integrated programmer permits fast and easy programming of setpoint temperature sequences. This temperature sequence is called program. A program is composed of individual sections (Step). The sections are defined by duration (t) or Gradient (°/t) and target temperature. The target temperature is the setpoint (SP), which is achieved at the end of a section. The programmer calculates the temperature ramp from the difference in time and temperature.

Create programs


A flashing segment indicates that a number needs to be entered. Under the submenu "Edit", enter a program number. Six programs may be stored (numbers 0 to 5).

When the program is running, only sections with complete information for the target temperature and time period are considered.

It makes sense to skip section numbers in the profile in order to use them later for corrections to the profile.

Important:

If a time of 00:00 is set for a profile, the profile is continued with the next section only after the setpoint temperature (±0.2 °C) is reached.

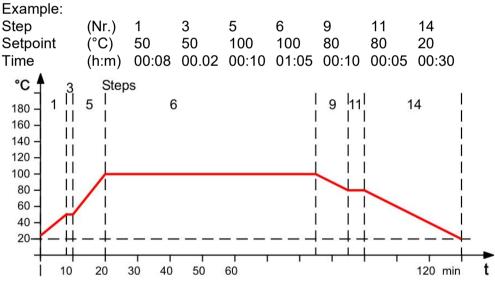

	Editing		
	PrgNr:	Set the program you wish to create. 0 5	
MENU V	SP :	Setpoint	
1 MAIN MENU ♣Start Program →Manage Programs ₽Pump OK ↓	TI : GRD: ST:	Time [hh:mm] or Gradient [°/t] = [°/min] Change type of program in MENU >configuration<, menu item >Programmtyp/program type< . Step 059	

(i) Press the	5	key if a	parameter is to	be retained.	(prior to OK)
---------------	---	----------	-----------------	--------------	---------------

	•		
Level 2 2 MANAGE PROGRAMS →Edit Pr9Nr. ▼Delete Pr9Nr. Print Pr9Nr.		Example: Program (PrgNr.) 2 is to have a target temperature with a set point (SW) of 100.00 °C in section t (ABS) 3. Duration (ZI) 1 hour.	
ок 🖡			
2 MANAGE PROGRAMS Edit \$ Pr9Nr.2) <u>/</u>	 Select the program number. Set the desired >PrgNr.< by pressing and οκ. 	
Level 3		For the selected section >ST< a setpoint >SP< and the duration >TI< or the gradient >GRD< are entered in the right half of the input window.	
3 EDITOR >Pr9Nr.2 SP :C 100.00C TI :T 1.00T ST : 2 ♣ 3		Input window for >ST<, >SP< and >TI< / >GRD< 1. Set the desired Step >ST < by pressing and οκ.	
or 3 EDITOR >Pr9Nr.2 SP :CC GRD:%% ST :		 Set set point >SP< by using the numeric keypad and confirm by pressing OK Set the time >TI< or a gradient >GRD< using the numeric keypad and confirm by pressing OK 	
(m)		Image by confirming time or GRD the values on the left side of the input window are moved.	
3 EDITOR >Pr9Nr.2 SP : 100.00CC TI : 1.00TT ST : 3 \$ 4	Ж	 The values for the next section are entered in the right half of the input window. See above 1. / 2. / 3. 	

Delete

It is possible to delete an entire program or consecutive sections

Press the 🕥 key if a parameter is to be retained. (prior to OK)

Delete PrgNr:	Set the program you wish to delete	
	0 5	
Delete from Step X	to Step X - (0 59)	

4	М	F	N	u	
Ν.	141	-		U	

1 MAIN MENU AStart Program →Manage Programs ▼Pump

Ψ

₩Pump				
ок 🖌				
Level 2		Submenu >Delete<		
2 MANAGE PROGRAMS ▲Edit Pr9Nr. →Delete Pr9Nr. ▼Print Pr9Nr.		 Set menu item >Delete PrgNr.< by pressing and ок 		
2 MANAGE PROGRAMS		 Set desired PrgNr with 		
Delete \$ Pr9Nr.0	Ж	Example: PrgNr. 0		
Level 3		Input window: delete sections.		
3 DELETE >Pr9Nr. 0 →from Step 0 ★to Step 59 delete 59 delete 59 from Step 0 from Step 0 from Step 0 J DELETE >Pr9Nr. 0 from Step 0 J DELETE >Pr9Nr. 0 to Step 59		 Select menu items > → from Step< and >→ to Step< one after the other by pressing and oκ. Select the desired section for every menu item by using the numeric keypad and confirm by pressing oκ. Example: >from 0< >to 59 		

		 Select menu item >delete< by pressing and ok
3 DELETE >Pr9Nr. 0 delete?	Ж	 Confirm >Delete?< by pressing Οκ
3 DELETE >Pr9Nr. 0 ! delete !		The delete procedure is confirm by indication of the message >delete< on the LCD display.

Print



ок 🗸

Every program can be printed via the serial interface for control and documentation purposes.

Press the key if a parameter is to be retained. (prior to OK)

Level 2	Submenu >Print<		
2 MANAGE PROGRAMS ★Delete Pr9Nr. →Print Pr9Nr.	 Set menu item >Print PrgNr.< by pressing and Οκ 		
2 MANAGE PROGRAMS Print \$ PrgNr.4	 Set desired >PrgNr.< by pressing and οκ Example: 4 		
Level 3			
3 PRINT Pr9Nr.4 →Print	• Confirm - → > Print< by pressing ок .		
3 PRINT Pr9Nr.4 Print?	 Confirm >Print?< by pressing ок . 		
3 PRINT PrgNr.4 ! print ! 3 PRINT PrgNr.4 done !	I Printing and end of printing are confirmed by successive messages on the LCD display as shown on the left.		

Setting the pump pressure 9.3.

MENU

1 MAIN MENU ▲Manage Programs →Pump **↓**Confi9uration ок 🕁

Ψ

The capacity of the circulating pump is set by adjusting the motor speed

① Press the ¹ key if a parameter is to be retained. (prior to OK)

Level 2		Submenu >Pump<	
2 PUMP →Pumpe sta9e 1		 The set pump stage is displayed. Example: 1 Confirm by pressing ок . 	
2 PUMP →Pumpe stage		 Set required pump stage by pressing and οκ Example: 3 	

Example:

Effective



Adjustable pump capacity stage 1 ... 4 Illuminated display: 🍣 for pump pressure

- (i) Indication of the adjusted pump pressure stage in the –OFF- mode.
- () Indication of the effective pump pressure stage (rotational speed) after start.

To protect the pump motor, the rotational speed changes with the load applied, for example, depending on the viscosity of the bath fluid at different working temperatures.

Flow rate:	22	26 l/min			
Pump capacity stage		1	2	3	4
Pump pressure	[bar] 0.4	0.5	0.6	0.7
Suction pump	[bar] 0.2	0.26	0.33	0.4
Total capacity	[bar] 0.6	0.76	0.93	1.1
in a loop circuit					

9.4. Configuration		
	2 CONFIGURATION	4
MENU	→Remote Setpoint ext Autostart Off-Mode p ActVar co Time/Date Language er	off off •. off ontrol
1 MAIN MENU ←Pump →Configuration →Control	Unit Program Type Reset C-Fühler 	°C Grad
Level 2 2 CONFIGURATION →Remote off *SetPoint ext off oK	Parameter level	 Remote control via serial interface. Possible parameters: off / on Switch on and off remote control by pressing First set the >type< of interface (RS232 or RS485) in the MENU >INTERFACE SERIAL< and then switch >Remote< to >on<. Control display in the top line for Remote ^o ^o ^o ^o Connect SERIAL with PC.
2 CONFIGURATION Aremote off →Setpoint ext off ★Autostart off Kor	 2 CONFIGURATION Setpoint ext #off Switch over setpoint setting by pressing and OK 	External setpoint setting Possible parameters: off / Eprog / PT100 • Switch over setpoint setting by pressing and ok off – Setpoint setting with the keys or Setpoint setting via the analog socket "ext. Pt100 or analog socket >REG+EPROG<
2 CONFIGURATION ★Setpoint ext off →Autostart off ▼Off-Mode p. off oK	2 CONFIGURATION Autostart \$ off	Autostart Possible parameters: on / off • Switch on and off Autostart by pressing and ok

	Parameter level	
2 CONFIGURATION Autostart off →Off-Mode P. off ▼ActVar control K	2 CONFIGURATION Off-Mode	 Pump-Mode Possible parameters: P. on / P. off Switch on and off >Off-Mode by pressing and ок P.on continuous operation of circulating pump P.off circulating pump is linked to Start/Stop
2 CONFIGURATION ★Off-Mode P. off →ActVar control ▼Time/Date →OK	2 CONFIGURATION ActVar \$ control	 ActVar - actuating variable Possible parameters: control, Eprog, serial ③ Set desired parameters with and OK. ④ Programming of variables for the parameters > serial < or > EPROG < is only accepted, if the unit is in Start mode
2 CONFIGURATION ▲ActVar control →Time/Date ▼Lan9ua9e en9lish		Time, Date, Year see next page
2 CONFIGURATION [★] Time/Date →Lan9ua9e en9lish ▼Unit °C ok	2 KONFIGURATION Lan9ua9e ∉en9lish	Language Possible parameters: deutsch / english • Set desired parameters with and OK.
2 CONFIGURATION ▲Lan9ua9e en9lish →Unit °C ▼Pro9ram type time	2 KONFIGURATION Unit	Unit Possible parameters: ° С / °F • Set desired parameters with and ок.
2 CONFIGURATION [♠] Unit °C →Program type time ▼Reset	2 KONFIGURATION Program type 🜩 time	 Program type Possible parameters: time / grad Set desired parameters with and OK .

Level 2	Parameter level	
2 CONFIGURATION [♠] Program type time →Reset oK	2 CONFIGURATION Reset? 2 CONFIGURATION Reset: Running	 Reset Return to factory settings by pressing ok >Reset?< returns all set values to the factory setting except for date and time. A RESET can be effected only in the – OFF- mode. During the message >Running < all parameters are reset to factory settings.
Level 2	Level 3	Parameter / Value
2 CONFIGURATION *ActVar control Jime/Date *Lan9ua9e en9lish CK	3 SET TIME/DATE →hour:min 11:50 Day.Mon 07.07 OK 3 SET TIME/DATE hour:min 11:50 →Day.Mon 07.07 Year 2009 CURRENT: 12:05:16 OK 3 SET TIME/DATE Year 2009 CURRENT: 12:05:16 OK 3 SET TIME/DATE Year 2009 CURRENT: 12:05:16 OK	Possible parameters: Time >hour:min Date >Day:Mon Year • Succesively set time, date and year using the numeric keypad and confirm by pressing OK . 3 SET TIME/DATE hour:min 11:50



The internal real time clock allows starting a program any time. The clock is set to the local mean time (MEZ) at the factory.

- (1) If the unit is operated in a different time zone, the clock can be adjusted in this menu.
- (i) Change summer/winter time in this menu

9.4.1. Remote control via the serial interface

Factory setting: off

,	0					
RS232	:	25.00°C				
IntAct	:	27.00°C				
Power	:	0%				
Control	:	intern				
RS485	:	25.00°C				
IntAct	:	27.00°C				
Power	:	0 %				
Control	:	intern				
0						

Setpoint is set via the serial RS232/RS485 interface through a PC or superordinated data system. In the header of the VFD COMFORT-DISPLAY, an "**R**" illuminates. It indicates that remote control mode is active.

← The selected setting is shown on the LCD DIALOG-DISPLAY.

o o	IMPORTANT: additional measures for remote control
SERIAL	(i) RS232 or RS485:
	First set >type< of interface in the MAIN MENU > Interface serial < the switch >Remote< to >on<.
	① Connect the circulator to the PC using an interface cable.
	Ocheck the interface parameters of both interfaces (circulator and PC) and make sure they match.
	(See 12.1. Setup for remote control page 82)

9.4.2. Keypad control or setpoint setting via the analog input

Factory setting: off

2	(:0	Ν	F	Ι	G	U	R	Α	Т	Ι	0	Ν			
-														 		
9	56	et	P	o	i	n	t		e	×	t			\$ 0	f	f
-														 		

The selected mode is indicated on standard display.

Setpoint	1:	152.00°C
IntAct	:	50.64°C
Power	:	100 %
Control	:	intern
Pt100	:	85.00°C
IntAct	:	83.00°C
Power	:	60 %
Control	:	external
Epro9		50.00°C
IntAct	1	27.40°C
Power	:	80 %
Control	:	internal

2 CONFIGURATION

→Remote		off	
Setpoint	ext	Epro9	
Autostart		off	
Off-Mode	P.	off	
ActVar		Epro9	
Time/Date			

In addition to the serial interface via remote control the circulator offers the possibility to adjust the setpoint via analog interface >ext. Pt100< or >REG+E-PROG<.

Possible parameters:

- **off** Setpoint is set via the keypad or via the integrated programmer. (factory setting
- **Pt100 -** Setpoint setting via the analog socket "ext. Pt100" using an external temperature sensor or an appropriate voltage/current source.
- **Eprog** Can only be adjusted when an electronic module with analog connections is used (option). Setpoint setting via the analog interface REG+E-PROG connection with an external voltage or current source or a programmer.

Important:

- Connect the external voltage or current source or a programmer to the circulator via the socket REG+E-PROG (see page 71).
 In the menu >MENU Inputs/Outputs< set the parameter >Erog-INPUT< and the input variables >Eprog-signal< (see page 71).
- The E-Prog input can only be used either under menu item >Setpoint ext < or under menu item >ActVar<</p>

9.4.3. Autostart

2 CONFIGURATION	
- Setpoint ext	off
→Autostart	off
♥Off-Mode P.	off

Possible parameters:

on - Autostart on

off - Autostart off (Factory setting)

Note:

The circulator has been configured and delivered by JULABO according to N.A.M.U.R. recommendations. This means for the start mode that the unit must enter a safe operating state after a power failure (non-automatic start mode). This safe operating state is indicated by "**OFF**" or "**R OFF**" on the VFD-COMFORT-DISPLAY. A complete shutdown of the main functional elements such as the heater and circulating pump is effected simultaneously.

The values set on the circulator remain stored, and the unit is returned to operation by pressing the start/stop key (in manual control mode). In remote control mode, the values need to be resent by the PC via the interface.

Should such a safety standard not be required, the AUTOSTART function (automatic start mode) may be activated, thus allowing the unit to be started directly by pressing the mains power switch or using a timer.

The AUTOSTART function can be used only if the setpoint is set via >keyboard<, >eprog<, or >ext. Pt100<.</p>



Warning

For supervised or unsupervised operation with the "AUTOSTART" function avoid any hazardous situation to persons or property

Take care to fully observe the safety and warning functions of the circulator.

9.4.4. Off-Mode

2 CONFIGURAT	[ON	
* Autostart		off
→Off-Mode	Ρ.	off
₩ActVar	cont	trol

OFF-Mode

Normally the circulating pump is switched via the start/stop signal. However, if circulation should be maintained even in the -off- condition, the parameter >**pump on<** needs to be set. Possible parameters:

- P.on Circulating pump in continuous operation
- P.off Circulating pump linked to Start/Stop (factory setting)

9.4.5. ActVar - actuating variable

2 CONFIGURAT	TON
daa sarsariii aa saatsariiiii	
# Off-Mode	P. off
→ActVar	control
∵ Time∕Date	

control

serial

Eprog

2 CONFIGURATION

2 CONFIGURATION

2 CONFIGURATION

ActVar

ActVar

ActUar

The variable (*ACTuating VARiable*) corresponds to the extent to which the heater or cooling unit of the circulator is controlled. Heat or cold is applied to the bath according to this variable. If this happens with the control electronics of the circulator, called > control < in this particular case, the bath temperature is exactly heated and maintained constant at the adjusted setpoint.

Programming of variables for the parameters > SERIAL < or > EPROG < is only accepted, if the unit is in Start mode.</p>

Possible parameters:

- **control** The internal control electronics of the circulator controls the heater and the connected cooling unit. Self-tuning is possible.. (factory setting)
- **serial** – The heater or the connected cooling unit receives the control signal via the serial interface. Self-tuning is not possible.
- **Eprog** The heater or the connected cooling unit receives the control signal via the E-Prog input. Self-tuning is not possible.
 - Setting requires electronic module.

Important:

Set the parameter >Eprog< and the input variables also in the menu >Inputs/Outputs< (see page 71).

2 CONFIGURATION

∙Remote		off
Setpoint	$e \times t$	Epro9
Autostart		off
Off-Mode	P.	off
ActVar		Epro9
Time/Date		

Note:

The E-Prog input can only be used either under menu item >Setpoint ext< or under menu item > ActVar <.



Warning:

The working temperature range of the circulator is determined during configuration. If set to >Control<, this range cannot be exceeded.

If set to > Serial < and > Eprog <, heat or cold is applied to the bath without control. The permissible maximum temperature can be exceeded. The user has to take adequate precautions for temperature control.

Materials, such as gaskets or insulations for example, may be damaged or destroyed, if the permissible maximum temperature is exceeded.

The safety and warning functions >

9.4.6. Setting of clock and date



2 CONFIGURAT	ION	
≜ ActVar	control	
→Time/Date		
▼Lan9ua9e	en9lish	

9.4.7. Language

The internal real time clock allows starting a program any time. The clock is set to the local mean time (MEZ) at the factory.

- (i) If the unit is operated in a different time zone, the clock can be adjusted in this menu.
- (i) Change summer/winter time in this menu

There are two options for the language of the LCD DIALOG-DISPLAY: German and English.

2 CONFIGURAT	ION
* Time/Date	
→Lan9ua9e	en9lish
‴Unit	°C

Possible parameters: deutsch / english

9.4.8. Unit

2 CONFIGURATI	ON.
*Lan9ua9e	english
→Unit	°C
♥Pro9ram typ	Zeit

Temperature values can be displayed in °C or °F. With this change, all adjusted temperature values are converted and displayed in the new unit. <u>Possible parameters:</u> °C (factory setting) °F

9.4.9. Program type

2 CONFIGURATION	
- Unit	°C
→Pro9ram typ	time
▼Reset	

A program can be created with the parameters time or gradient. <u>Possible parameters:</u> **time** (factory setting) **Grd** (Gradient) [°/t] = [°/min]

9.4.10. Reset

2 CONFIGURATION	
♣Pro9ram typ	time
→Reset	

A Reset will return all values to factory setting except for date and time.

(i) A Reset can be effected in the >OFF< mode only.

9.5. Control

The circulator is designed for internal and external temperature control. This can be switched in this submenu. Only the set of parameters corresponding to the particular setting is displayed.

	Control internal		trol external	
	<pre> →ControlType ▼Selftunin9 Dynamic Xp Tn Tv </pre>	always ▼Se	n 720s 9 55s	
MENU 1 MAIN MENU Configuration Control Interface serial OK	 For external temper external sensor to the Press the key 	ne socket at the r	ear of the circulator.	

Level 2	Parameter-Level	
2 CONTROL →ControlType inter ▼Selftunin9 alway:		 Switch over control type by pressing and οκ
);;;;;;;;;;;;;;;;;;;;;;;;;;;;;;;;;;;;;	Possible parameters: intern / extern
	Or 2 CONTROL ControlType textern	 The control type can be adjusted in the -OFF- mode only.
2 CONTROL ★ControlType intern →Selftunin9 always ★Dynamic aperic	2 CONTROL Selftuning always	 The parameter flashes, switch by pressing and ok Possible parameters: off / once / always.
Control Type internal	Parameter-Level	
2 CONTROL		① The parameter flashes, switch by pressing

Control Type internal	Parameter-Level	
2 CONTROL ▲Selftunin9 always →Dynamic aperio ▼XP 1.0K	2 CONTROL Dynamic aperio	 The parameter flashes, switch by pressing and οκ Possible parameters: aperio / normal This parameter affects the temperature sequence in case of internal control.

	Parameter-Level	
2 CONTROL Denamik aperio →XP 1.0K Th 100s	2 CONTROL Xp 1.0K	 The value flashes. Set by using the numeric keypad and confirm by pressing ок. Example: 1 т 0 ок Setting range: 0.1 99.9 К
2 CONTROL [▲] XP →Tn 1.0K →Tn 100s ▼Tv 5s OK	2 CONTROL 	 The value flashes. Set by using the numeric keypad and confirm by pressing ок. Setting range: 3 9999 s
2 CONTROL ▲Tn 100s →Tv 5s 	2 CONTROL 	 The value flashes. Set by using the numeric keypad and confirm by pressing Οκ.
Control Type external	Parameter-Level	
2 CONTROL ★Selftunin9 always →CoSpeed 0.0 ▼XP 0.7K	2 CONTROL CoSpeed 0.0	 The value flashes. Set by using the numeric keypad and confirm by pressing ок. Example: 0 т 0 ок
		Setting range: 0.0 5.0
2 CONTROL ★CoSpeed 0.0 →XP 0.7K ▼Tn 720s	2 CONTROL XP 0.7K	 The value flashes. Set by using the numeric keypad and confirm by pressing ок. Setting range: 0.1 99.9 К
2 CONTROL ▲XP 0.7K →Tn 720s ▼T∨ 55s	2 CONTROL Tn 720s	 The value flashes. Set by using the numeric keypad and confirm by pressing ок. Setting range: 3 9999 s
2 CONTROL ▲Tn 720s →Tv 55s ▼XPu 2.9K	2 CONTROL Tv 55s	 The value flashes. Set by using the numeric keypad and confirm by pressing ок. Setting range: 0 999 s
2 CONTROL ▲Tv 55s →Xpu 2.9K 	2 CONTROL Xpu 2.9K	 The value flashes. Set by using the numeric keypad and confirm by pressing ок. Setting range: 0.1 99.9 К

9.5.1. Control internal / external



Pt100

(i) The control type can be adjusted in the **-OFF-** mode only.

Possible parameters:

internal temperature control

extern external temperature control with external Pt100 sensor

IMPORTANT:

Additional measures for external temperature control.

- (i) Connect a Pt100 sensor to the socket on the rear of the circulator.
- Sensor calibration of the external Pt100 sensor is performed in the >ATCalibration< menu.</p>
- Suggested adjustments for external temperature control: Upper band / Lower band and IntMax / IntMin see Section >Limits<.</p>
- External temperature control does not allow the setpoint to be set via the "ext. Pt100" socket.



Notice:

Place the external sensor into the temperature-controlled medium and securely fix the sensor.

Accessory: Pt100 external sensor

Order No.	Description	Material	Cable
8981003	200x6 mm Ø,	stainless steel	1.5 m
8981005	200x6 mm Ø,	glass	1.5 m
8981006	20x2 mm Ø,	stainless steel	1.5 m
8981010	300x6 mm Ø,	stainless steel	1.5 m
8981015	300x6 mm Ø,	stainless steel / PTFE coated	3 m
8981013	600x6 mm Ø,	stainless steel / PTFE coated	3 m
8981016	900x6 mm Ø,	stainless steel / PTFE coated	3 m
8981014	1200x6 mm Ø,	stainless steel / PTFE coated	3 m
8981103	Extension cable for	Pt100 sensor	3.5 m
8981020	M+R in-line Pt100	sensor	

Pt100

M+R

R

m Ab

The M+R in-line Pt100 sensor is a flow sensor and can be installed loop circuit

9.5.2. Selftuning

When performing a selftuning for the controlled system (temperature application system), the control parameters Xp, Tn and Tv are automatically determined and stored.

Possible parameters:

off - <u>no</u>selftuning

The control parameters ascertained during the last identification are used for control purposes.

once - single selftuning (factory setting)

The instrument performs a single selftuning of the controlled system after each start with the ok key or after receiving a start command via the interface.


always - continual selftuning

The instrument performs a selftuning of the controlled system whenever a new setpoint is to be reached.

Use this setting only when the temperature application system changes permanently.

9.5.3. Dynamik intern

This parameter affects the temperature sequence only in case of internal control.

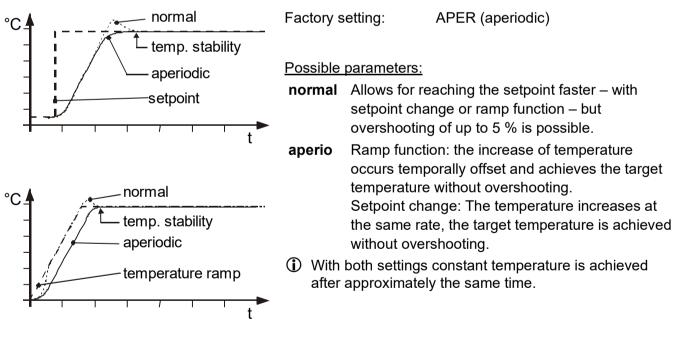


9.5.4. CoSpeed - extern

CoSpeed:

This parameter affects the temperature pattern only in case of **external** control.

Possible parameters: 0 ... 5

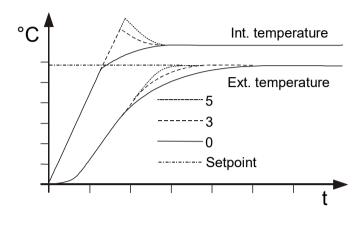

During selftuning, the control parameters Xp, Tn, and Tv of a controlled system are automatically determined and stored. The time required for tuning may vary depending on the controlled system. This controller design allows protection of sensitive objects requiring temperature control. As soon as a co-speed factor is set, it is considered in calculating the control parameters. As shown in the diagram, tuning times are shorter the higher the co-speed factor is, but overshoot can occur in the internal system.


9.5.5. Control parameters – internal/external

Control parameters - internal/external

The control parameters preset at the factory are in most cases adequate to achieve an optimal temperature pattern in the samples requiring temperature control.

Each parameter may be manually set via the keypad if necessary to allow optimal control performance.

Setting range:	Proportional range >Xp<
internal / external 0.1 99.9 K	The proportional range is the temperature range below the setpoint temperature value in which the control circuit reduces the heating power from 100% to 0%.
Setting range:	<u>Resetting time >Tn<</u> (Integral component)
internal / external 39999 s	Compensation of the remaining control deviation due to proportional control. An insufficient resetting time may cause instabilities to occur. An excessive resetting time will unnecessarily prolong compensation of the control difference.
Setting range:	Lead time >Tv< (Differential component)
internal / external 0 999 s	The differential component reduces the control settling time. An insufficient lead time will prolong the time required to compensate for disturbance effects and cause high overshoot during run-up. An excessive lead time could cause instabilities (oscillations) to occur.
Setting range: 0.1 99.9 K	<u>Proportional range >Xpu<</u> The proportional range Xpu of the cascaded controller is only needed for external control.

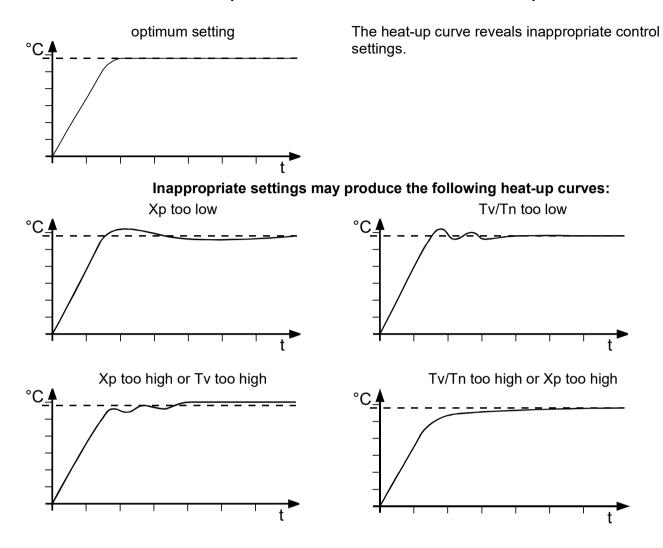

Optimization instructions for the PID control parameters:

9.6. Serial Interface

Factory settings: 4800 Baud even hardware handshake For communication between the circulator and a PC or a superordinated process system, the interface parameters of both units must be identical. The adjustment is usually carried out one time only in the >Interface serial< menu.

RS485

1 MAIN MENU •Control	
→Interface serial ▼ATCalibration	
ок 🗸	

2 INTERFACE	SERIAL
→Type ▼Baudrate	RS232 4800
Parity	even
Handshake	hardw

2 INTERFACE	SERIAL
÷Туре	RS485
▼Baudrate	4800
Parity	even
Handshake	hardw
Adress	0

Level 2	Parameter-level	Press the key if a parameter is to be retained. (prior to OK)
2 INTERFACE SERIAL →Type RS232 *Baudrate 4800 OK →	2 INTERFACE SERIAL Type	 The parameter flashes, switch by pressing and ок Possible parameters: RS232 / RS485 Address RS485: 0 to 127
2 INTERFACE SERIAL [♠] Type RS232 →Baudrate 4800 ▼ Parity even OK →	2 INTERFACE SERIAL Baudrate \$ 4800	 The parameter flashes, switch by pressing and ok Possible parameters: 4800 / 9600 / 19200 / 38400
2 INTERFACE SERIAL ★Baudrate 4800 →Parity even ▼Handshake hardw OK →	2 INTERFACE SERIAL Parity \$ even	 The parameter flashes, switch by pressing and ok Possible parameters: no/even/odd even: Data bits = 7; Stop bits = 1 odd: Data bits = 7; Stop bits = 1 no: Data bits = 8; Stop bits = 1
2 INTERFACE SERIAL [★] Parity even →Handshake hardw 	2 SERIELLES INTERF. Handshake 🖨 hardw	 The parameter flashes, switch by pressing and ок Possible parameters: hardw / softw Xon/Xoff-protocol (Software handshake) Protocol RTS/CTS (Hardware handshake)

Example:


ATC Absolute Temperature Calibration, 3-point calibration 9.7.

ATC is used to compensate for a temperature difference between the circulator and a defined measuring point in the bath tank that may develop for physical reasons.

internal ATC

internal			external		
2 ATCalibra	ət	ion	2 ATCalibra	at	ion
→Sensor		intern	⇒Sensor		extern
▼Status		no	▼Status		no
Туре		3-point	Туре		3-point
TmpValue1	:	××.××°C	TmpValue1	:	××.××°C
CalValue1	:	××.××°C	CalValue1	:	××.××°C
TmpValue2	:	××.××°C	TmpValue2	:	××.××°C
CalValue2	:	××.××°C	CalValue2	:	××.××°C
TmpValue3	:	××.××°C	TmpValue3	:	××.××°C
CalValue3	:	×ו××°C	CalValue3	i	××.××°C

Principle: For ATC calibration, in steady state the bath temperature at the location of the temperature sensor (CT) is determined at the respective adjusted Μ

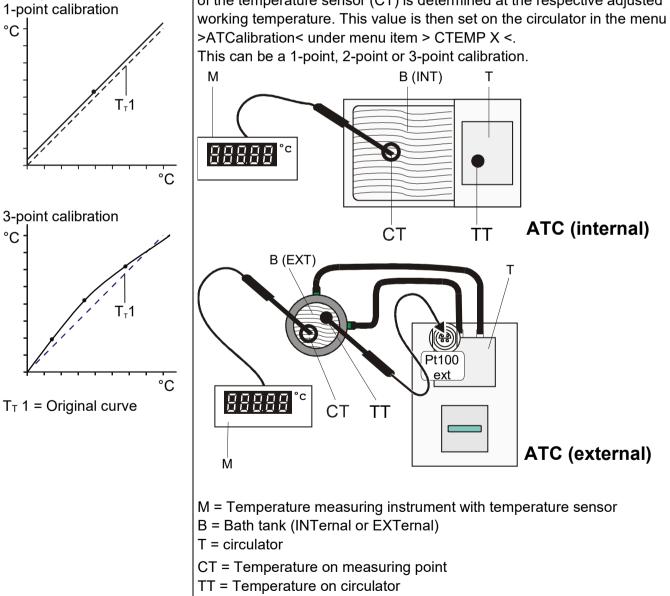

MENU 1 MAIN MENU ▲Interface serial →ATCalibration ▼Limits OK	 Press the <a>key if a 	parameter is to be retained. (prior to OK)
Level 2	Parameter-Level	
2 ATCalibration →Sensor intern ▼Status no OK	2 ATCalibration Sensor intern	 The parameter flashes, switch by pressing and
2 ATCalibration * Sensor intern	2 ATCalibration	The parameter flashes, switch by pressing
→Status no ▼Type 3-point OK	or 2 ATCalibration 	 and ок >no< Carry out an ATC calibration)yes< return to standard operation after calibration.
2 ATCalibration ▲Status no →Type 3-point ▼TmpValue1 : xx.xx°C OK	2 ATCalibration Type 1-Point	 The parameter flashes, switch by pressing and οκ A >1-point<, >2-point< or >3-point<
	Type 2-point 2 ATCalibration Type 3-point	calibration can be carried out.
2 ATCalibration *Type 3-Point →TmpValue1 : 80.00°C ▼CalValue1 : 79.70°C OK	2 ATCalibration TmpValue1 : xx.xx°C	 The value >TmpValue< is only indicated In addition the measured temperature value >CalValue X< is saved during the next step.
2 ATCalibration [▲] TmpValue1 : xx.xx [®] C →CalValue1 : xx.xx [®] C ▼TmpValue2 : xx.xx [®] C OK	2 ATCalibration CalValue1 : xx.xx°C	 The value flashes. Set by using the numeric keypad and confirm by pressing οκ
	 If only a 1-point calibration not indicated anymore 	on is carried out, the following menu items are

2 ATCalibration *CalValue1 : xx.xx°C →TmpValue2 : xx.xx°C *CalValue2 : xx.xx°C OK	2 ATCalibration TmpValue2 : xx.xx°C	 The value >TmpValue< is only indicated In addition the measured temperature value >CalValue X< is saved during the next step.
2 ATCalibration [★] TmpValue2 : xx.xx°C →CalValue2 : xx.xx°C ▼TmpValue3 : xx.xx°C OK	2 ATCalibration CalValue2 : xx.xx*C	 The value flashes. Set by using the numeric keypad and confirm by pressing οκ
	If only a 2-point calibratio not indicated anymore	n is carried out, the following menu items are
2 ATCalibration *CalValue2 : xx.xx°C →TmpValue3 : xx.xx°C *CalValue3 : xx.xx°C OK	2 ATCalibration TmpValue3 : xx.xx°C 	 The value >TmpValue< is only indicated. In addition the measured temperature value >CalValue X< is saved during the next step
2 ATCalibration [★] TmpValue3 : xx.xx°C →CalValue3 : xx.xx°C OK	2 ATCalibration CalValue3 : xx.xx°C	 The value flashes. Set by using the numeric keypad and confirm by pressing οκ

9.7.1. ATC Fühler - intern / extern

2 ATCalibra	tion
→Sensor	intern
▼Status	no

In the first submenu the ATC function is set for the >intern< internal or the >extern< external temperature sensor.

Calibration can be carried out for the internal temperature sensor and for the external temperature sensor connected to the socket "ext. Pt100". The circulator is able to save both parameter sets. However only the one which has been set under menu item > sensor < is displayed.

9.7.2. ATC Status - yes / no

2 ATCali	bration	
≜ Sensor	intern	
⇒Status	no	
₩Туре	3-point	

In the second submenu the ATC function for the temperature sensor selected above is activated >yes< or deactivated >no<.

>yes< (factory setting) The controller of the circulator uses the original curve of the temperature sensor or the new curve measured during the ATC calibration.


Important: Set to >NO< during the calibration process

>no< An ATC calibration is to be carried out. Important: Set to >yes< after calibration.</p>

(i) In the > Status < >yes< the ATC calibration always affects the current working temperature; also the one set via interface.

9.7.3. Type 1 -/ 2 -/ 3 point

2 ATCalibra	tion
Туре	1-point
2 ATCalibra	tion
Туре 	2-point
2 ATCalibra	tion
Туре	3-point

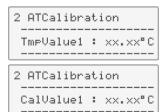

A >1-point<, >2-point< or >3-point< calibration can be carried out.

First geometrically define the location for calibration (measuring point CT), then determine the temperature values of the calibration points. The type of calibrations also determines the number of the following pairs of values indicated on the LCD DIALOG-DISPLAY.

Pairs of values:

TmpValue X: Circulator temperature 1 or 2 or 3 (actual value TT) The actual temperature of the bath is simultaneously saved with the "calibration value" >CalValue< and can be indicated for control purposes (value does not flash).

CalValue X: Calibration temperature 1 or 2 or 3 (actual value CT) The "calibration value" is determined with a temperature measuring device and saved under menu item >CalValue<. (value flashes (1))

9.7.4. Example: 3-point calibration for internal control

In the temperature range from 80 $^{\circ}$ C to 160 $^{\circ}$ C the calibration curve of the temperature sensor (TT) is to be adjusted to the actual temperatures at measuring point (CT).

2 CONTROL →ControlType intern *Selftuning always OK 80.00 °C 120.00 °C 120.00 °C 160.00 °C SetPoint1: \$80.00°C IntAct : 24.64°C Power : 0 % Control : internal	 2. Set working temperature set See standard display line 1 Example 80.00 °C Press the keys 8 and The bath is heated up. 	set only in the –OFF- mode.	
Pt100 Pt100 79.70 °C ct	3. Reading of temperature measuring device Read the value of measuring point CT on the device and enter under menu item > CalValue X< by using the keypad. 		
MENU	4. Calibration		
1 MAIN MENU ▲Interface serial →ATCalibration ▼Limits OK	 Press the key if a parameter is to be retained. (prior to OK) Setting is required only for the first calibration point. 		
Level 2	Parameter-Level		
2 ATCalibration →Sensor intern ▼Status no	2 ATCalibration Sensor intern	 Set sensor: >intern<. The parameter flashes, switch by pressing and οκ 	
2 ATCalibration ▲Sensor intern →Status no ▼Type 3-point OK	2 ATCalibration Status no	 Set to >no The parameter flashes, switch by pressing oκ 	
2 ATCalibration ▲Status no →Type 3-point ▼TmpValue1 : xx.xx°C OK	2 ATCalibration Type 3-point	 A >3-point< calibration is carried out. The parameter flashes, switch by pressing and οκ 	
2 ATCalibration AType 3-Point →TmpValue1 : 80.00°C ▼CalValue1 : 79.70°C OK	2 ATCalibration TmeValue1 : 80.00°C	 The value >TmpValue< is only indicated. In addition the measured temperature value >CalValue X< is saved during the next step. 	

2 ATCalibration [★] TmpValue1 : 80.00°C →CalValue1 : 79.70°C ▼TmpValue2 :120.00°C OK	2 ATCalibration CalValue1 : 79.70°C	 The value flashes. Set by using the numeric keypad and confirm by pressing ок
	Return to 2. Set working temp	erature value: 120.00 °C
2 ATCalibration ▲CalValue1 : 79.70°C →TmpValue2 :120.00°C ▼CalValue2 :119.50°C OK	2 ATCalibration TmpValue2 :120.00°C	 The value >TmpValue< is only indicated. In addition the measured temperature value >CalValue X< is saved during the next step.
2 ATCalibration ★TmpValue2 :120.00°C →CalValue2 :119.50°C ▼TmpValue3 :160.00°C OK	2 ATCalibration CalValue2 :119.50°C	 The value flashes. Set by using the numeric keypad and confirm by pressing OK .
	Return to 2. Set working temp	erature value: 160.00 °C
2 ATCalibration *CalValue2 :119.50°C →TmpValue3 :160.00°C ▼CalValue3 :159.30°C	2 ATCalibration TmeValue3 :160.00°C	 The value >TmpValue< is only indicated. In addition the measured temperature value >CalValue X< is saved during the next step.
2 ATCalibration [★] TmpValue3 :160.00°C →CalValue3 :159.30°C OK	2 ATCalibration CalValue3 :159.30°C	 The value flashes. Set by using the numeric keypad and confirm by pressing οκ
	5. Return to standard operat	ion
2 ATCalibration ▲Sensor intern →Status no ▼Type 3-point OK	2 ATCalibration Status yes	Set >YES< after calibration. (Standard operation) • The parameter flashes, switch by pressing and ок

9.8. Limits

The >Limits< menu allows the minimum and maximum values to be set for all important setting ranges and capacity variables.

Control internal		Control external
	2 LIMITS	
	→SetMax 300. →SetMin -94. Heatin9Max Coolin9Max	 .00°C →SetMax 300.00°C .99°C ▼SetMin -94.99°C 100% Heatin9Max 100%
MENU		
1 MAIN MENU ATCalibration →Limits ▼Inputs/Outputs OK	 Press the Skey if 	a parameter is to be retained. (prior to OK).
Level 2	Parameter-level	
2 LIMITS →SetMax 300.00°C ▼SetMin -94.99°C OK	2 LIMITS SetMax 300.00°C	 The parameter flashes. Set by using the numeric keypad and confirm by pressing OK
2 LIMITS [▲] SetMax 300.00°C →SetMin -94.99°C ▼Heatin9Max 100% OK	2 LIMITS SetMin -94.99°C	 The parameter flashes. Set by using the numeric keypad and confirm by pressing οκ
2 LIMITS [♠] SetMin -94.99°C →Heatin9Max 100% ▼Coolin9Max 0% OK	2 LIMITS HeatingMax 100%	 The parameter flashes. Set by using the numeric keypad and confirm by pressing οκ
2 LIMITS [♠] Heatin9Max 100% →Coolin9Max 100% ▼IntMax 300.00°C OK	2 LIMITS CoolingMax 0%	 The parameter flashes. Set by using the numeric keypad and confirm by pressing οκ
2 LIMITS ★Coolin9Max 0% →IntMax 300.00°C ▼IntMin -94.99°C OK	2 LIMITS IntMax 300.00°C	 The parameter flashes. Set by using the numeric keypad and confirm by pressing ок

2 LIMITS ▲IntMax 300.00°C →IntMin -94.99°C ▼Upper band 200K	2 LIMITS IntMin -94.99°C	•	The parameter flashes. Set by using the numeric keypad and confirm by pressing οκ .
2 LIMITS ▲IntMin -94.99°C →Upper band 200K ▼Lower band 200K	2 LIMITS UPPer band 200K	•	The parameter flashes. Set by using the numeric keypad and confirm by pressing OK .
2 LIMITS ▲Upper band 200K →Lower band 200K	2 LIMITS Lower band 200K	•	The parameter flashes. Set by using the numeric keypad and confirm by pressing or .

9.8.1. Limits for internal control and external control

2 LIMITS	
SetMax	300.00°C
2 LIMITS	
SetMin	-94.99°C

Set Max / Min– Maximum and minimum setpoint Restriction of the adjustable temperature range.

The limitation of the operating temperature range effects the temperature

setting in the menu with the key

Only setting of working temperatures which lie within the determined limits is possible

Existing settings for Setpoint 1, -2, -3, as well as those for >OverTemp< and > SubTemp <, are automatically deferred into the limit range.

Setting range: -94,90 °C ... +300,0 °C

2 LIMITS
Heatin9Max 100%
2 IMITS
Coolin9Max 0%

Set maximum heating / cooling

The heating and cooling capacity of the unit are adjustable. 100 % corresponds to the technical specification of the equipment.

Setting range:

Heating Max –	0 to 100 % in 1 % steps
Cooling Max –	0 to 100 % in 1 % steps

9.8.2. Limits for external control

2 LIMITS	
IntMax	300.00°C
2 LIMITS	
IntMin	-94.99°C



2 LIMITS

2 LIMITS

Upper band

Lower band

200K

200K

INTERN MAX / MIN Restriction for the temperature range of the internal bath.

Setting range: -94,9 °C ... +300,0 °C

The limits IntMax and IntMin are only active in external control. IntMax and IntMin determine fixed limits for the temperature within the internal bath. The temperature controller cannot exceed these limits even if it would be necessary for achieving the temperature in an external system. Therefore it is possible that the external setpoint cannot be achieved. Sense of limit setting:

Sense of limit setting:

- \blacksquare Protects the bath fluid from overheating.
- ✓ Prevents an undesired alarm shutdown by the excess temperature protection >ALARM CODE 14<.</p>

Set the value of > Int Max at least 5 °C below the value of >SafeTemp<.


- Protects the pump motor from high viscosity of the bath fluid at low temperatures.
- For refrigerated circulators. Freezing protection when using water as bath fluid.

Upper band: and Lower band:

The band limitation is active during external control. Varied, practiceoriented setting are feasible for heat-up and cool-down phases.

Setting range: 0 °C ... 200 °C


Upper band: and **Lower band** allow for the limitation of the difference between the temperatures in the internal bath and the external system to any maximum value for the heat-up and cool-down phase. During the heat-up phase this difference value is always added to the actual external temperature. During the cool-down phase the difference value is subtracted.

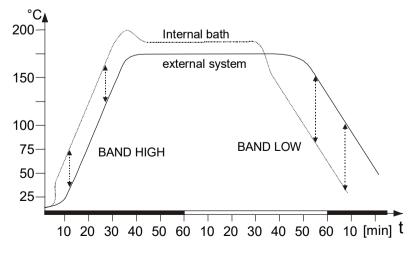

Sense of a band limitation:

- Protection of objects and samples by gentle temperature control
- Protection of e.g. glass reactors from thermal shock.

9.9. Analog inputs/outputs

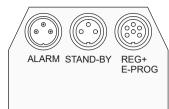



C-Pulse

C-Pulse ActInt

| Setpoint

inactiv

StdBy

Power Setpoint

2 INPUTS/OUTPUTS

2 INPUTS/OUTPUTS

- --- --- ---- ----

SV-OUT

→SV-OUT

Chan3

EPROG Ext.StBy

AlarmOut |

<mark>∞Chan1</mark> Chan2

(i) In order to use the analog inputs and outputs, the circulator must be	•
equipped with the electronic module, which is available as option.	
Order No. 8 900 100 Electronic module	

This submenu enables the input and output values to be set for the programmer input and the temperature recorder outputs of the REG+E-PROG socket.

The >STAND-BY< input and the >ALARM< output are configurable.

← If the electronic module is not installed, this menu can be used only to switch the control of the solenoid valve (SV).

SV-Output	Control connector (14) 230 V max. 1.25 A		
Channel 1	voltage output for recorder (V)		
Channel 2	voltage output for recorder (V))		
Channel 3	current output for recorder (mA)		
EPROG	external programmer input		
External standby Standby input			
	(for external switch-off)		
Alarm output	Output for external alarm signal		

MENU ↓ 1 MAIN MENU ▲Limits →Inputs/Outputs 		d for the connection of an HST Booster Heater. a parameter is to be retained. (prior to OK)
Level 2	Parameter-Level	
2 INPUTS/OUTPUTS -SV-OUT C-Pulse *Chani ActInt	2 INPUTS/OUTPUTS SV-OUT & C-Pulse Or 2 INPUTS/OUTPUTS SV-OUT & Backflow	Control output (14) 230 V max. 1,25 A: Possible parameters: C-Pulse, Backflow, Refill • The parameter flashes, switch by pressing and ок
	or 2 INPUTS∕OUTPUTS SV-OUT ♥ Refill	

Level 2	Level 3	Parameter-level
2 INPUTS/OUTPUTS ★SV-OUT C-Pulse →Chan1 ActInt ▼Chan2 Power	3 Channel1 →OutPut ActInt ♥ 0U = -99.90°C 10U = 400.00°C 3 Channel1 →OutPut ActInt	Define the output value for channel 1 Possible parameters: ActInt, ActExt, Power, Setpoint • The parameter flashes, switch by pressing and 3 Channel1
	3 Channel1 *0U *0utput ActInt *0U *0U <t< td=""><td>OutPut * ActInt Select the scale for channel 1. • The value flashes. Set by using the numeric keypad and confirm by pressing OK 3 Channel1 0U = -99.90°C</td></t<>	OutPut * ActInt Select the scale for channel 1. • The value flashes. Set by using the numeric keypad and confirm by pressing OK 3 Channel1 0U = -99.90°C
	3 Channel1 ★0V = -99.90°C →10V = 400.00°C OK	 The value flashes. Set by using the numeric keypad and confirm by pressing ΟΚ . 3 Channel1 10V = 400.00°C
2 INPUTS/OUTPUTS Chan1 ActInt Chan2 Power Chan3 SetPoint Chan3 SetPoint	3 Channel2 →Output Power ▼0V = 0.00 % 10V = 100.00 % 3 Kanal2 →Output Power ▼0V = 0.00 % ▼K	Define the output value for channel 2. Possible parameters: ActInt, ActExt, Power, Setpoint • The parameter flashes, switch by pressing and OK 3 Channel2 Output Power
	3 Channel2 ★OutPut Power →0V = 0.00 % ▼10V = 100.00 % ▼ ▲	Select the scale for channel 2. The value flashes. Set by using the numeric keypad and confirm by pressing OK . Channel2 OU = 0.00 %
	3 Channel2 ★0V = 0.00 % →10V = 100.00 % 	The value flashes. Set by using the numeric keypad and confirm by pressing OK .

Level 2	Level 3	Parameter-level
2 INPUTS/OUTPUTS [★] Chan2 Power →Chan3 SetPoint ▼EPROG SetPoint ok	3 Channel3 →Output Setpoint ▼0mA = -99.90°C 20mA = 400.00°C Region 0-20mA 3 Channel3 →Output Setpoint	Define the output value for channel 3 Possible parameters: ActInt, ActExt, Power, Setpoint • The parameter flashes, switch by pressing and OK 3 Channel3
	▼0mA =OK	Output \$ SetPoint
	3 Channel3 ▲Outrut Setroint →0mA = -99.90°C ▼20mA = 400.00°C	 The value flashes. Set by using the numeric keypad and confirm by pressing OK 3 Channel3 OmA = -99.90°C
	3 Channel3 ★0mA = -99.90°C →20mA = 400.00°C ▼Region 0-20mA OK	 The value flashes. Set by using the numeric keypad and confirm by pressing OK. 3 Channel3 > 20mA = 400.00° C
	3 Channel3 ★20mA = 400.00°C →Region 0-20mA oK	Possible parameters: 0-20 mA, 4-20 mA • The parameter flashes, switch by pressing and ok 3 Channel3 Region \$ 0-20mA
2 INPUTS/OUTPUTS Chan3 Setpoint →EPROG Setpoint Ext.StBy inactiv	3 EPROG →Input Setpoint Signal Voltage L Value -99.90°C H Value 400.00°C Setpoint : -xx.xx°C 	Possible parameters: Setpoint, variable, flow rate, pressure • The parameter flashes, switch by pressing and OK 3 EPROG Input \$ SetPoint 3 EPROG Input \$ Variable 3 EPROG Input \$ Flowrate 3 EPROG Input \$ Pressure

Level 2	Level 3	Parameter-level
	3 EPROG Input Setpoint →Signal Voltage L Value -99.90°C OK	Possible parameters: voltage, current The parameter flashes, switch by pressing and 3 EPR0G Signal \$ Voltage Vitage Signal \$ Current
	3 EPROG Signal Volta9e →L Value -99.90°C H Value 400.00°C	 The value flashes. Set by using the numeric keypad and confirm by pressing OK 3 EPROG L Value -99.90° C Viv
	3 EPROG L Value -99.90°C →H Value 400.00°C Setpoint : -xx.xx°C OK	 The value flashes Set by using the numeric keypad and confirm by pressing ок. 3 EPROG H Value 400.00° С
	3 EPROG H Value 400.00°C →Setpoint : -xx.xx°C oK	The value >setpoint< is only indicated. 3 EPR0G Setpoint : -xx.xx*C
2 INPUTS/OUTPUTS EPROG Setpoint →Ext.StBy inactiv AlarmOut StdBy	(Parameter Level) 2 INPUTS/OUTPUTS →Ext.StBy ♥ activ 2 INPUTS/OUTPUTS →Ext.StBy ♥ inactiv	 Possible parameters: inactive, active The parameter flashes, switch by pressing and οκ
2 INPUTS∕OUTPUTS Ext.StBy inactiv →AlarmOut StdBy 	3 ALARM OUTPUT 	Possible parameters: Alarm, Stand By, Alarm+ Stand By • The parameter flashes, switch by pressing ok and 3 ALARM OUTPUT • Function $=$ Alarm ·

Level 2	Level 3	Parameter-level
Level 2	Level 3 3 ALARM OUTPUT Function StdBy →TYPE normal 	Parameter-level Possible parameters: normal, inverse The parameter flashes, switch by pressing and ok ALARM OUTPUT TYPE TYPE TYPE TYPE TYPE TYPE TYPE TYPE TYPE Type Typ

9.9.1. Control connector



Control connector (14):


- SV-Out >Backflow< or >C-Pulse< or >Refill<
 - The control output (14) can be adapted to solenoid valves for different tasks.
- >Backflow< In the OFF condition, the control output (14) is not powered.</p>

See - Backflow safety device- on page 21

- >C-Pulse< The control output (14) can be used to emit a cooling pulse. For the pulsing period, the output is powered. (only used with Heating Circulators)
- >**Refill**< Control of a Automatic Refill Device (Accessories)

9.9.2. Outputs of the connector - REG+E-PROG

Outputs of the connector

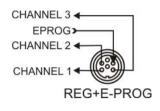

1. First define the desired output value for channels 1 to 3:

ActIntinternal actual temperature value (bath temperature)ActExtexternal actual temperature value (external sensor)Powerperiodic or intermittent heating or coolingSetpointactive setpoint temperature(setpoint 1, 2, 3/integr. programmer/external programmer)

2. Then select the scale for channels 1 to 3:

Voltage outputs channels 1 and 2


Assign the voltage values of 0 V to the lowest and 10 V to the highest necessary temperature or power rating required as an output value ($^{\circ}C/_{\%}$).

Current output channel 3

Examples:

Assign the current values 0 mA or 4 mA to the lowest and 20 mA to the highest temperature or power rating required as an output value ($^{\circ}C/\%$).

3. The current output (channel 3) offers 2 ranges for selection: 0 to 20 mA or 4 to 20 mA

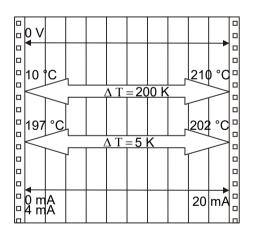

3 Char	3 Channel1 ≜Out⊧ut ActInt		
≜Outpu	ıt	ActInt	
÷ØŲ	=	10.00°C	
▼ 100	=	210.00°C	

lowest temperature value: 10 °C highest temperature value 210 °C Fig. shows 200 °C scaled to paper width slope: 50 mV/°C

lowest temperature value: 197 °C highest temperature value: 202 °C Fig. shows 5 °C scaled to paper width slope: 2000 mV/°C

9.9.3. Input of the connector - REG+E-PROG

E-PROG - Input

Menu >Configuration<

	ing congarate	•
	2 CONFIGURATIO	IN
	* Remote	off
	→Setpoint ext	Epro9
	▼Autostart	off
	2 CONFIGURATIO	IN
	≜ Off-Mode	P. off
	→ActVar	Epro9
2.	∵ Time∕Date	

Menu >Inputs/Outputs<

2 INPUTS/OU	JTPUTS
Chan3	Setpoint
→EPROG	Flowrate
Ext.StB9	inactiv

2 CONFIGURATION

Remote		off
Setpoint	ext	Epro9
Autostart		off
Off-Mode	P.	off
ActVar		Epro9
Time/Date		

Setting is necessary if

1. the Setpoint is to be set via an external voltage or current source or programmer

For this, in the menu >Configuration<, first set the menu item >Setpoint< to >Eprog<.

2. the heater variable should be controlled via an external control pulse.

For this, in the menu >Configuration<, set the menu item > ActVar.< to >Eprog<.

- 3. the signal of an external flow meter should be registered.
- (i) The E-Prog input can only be used either under menu item >Setpoint< or under menu item > Actvar <. If the input is occupied neither by >Setpoint< or > ActVar <, the signal of a flow meter can be connected.
- () First set menu >Configuration<, then menu >Inputs/Outputs<.
- Connect the external voltage or current source or programmer to the REG+E-PROG socket (12) of the circulator.

Selecting the signal:

The programmer (E-PROG) input of the circulator can be matched to the output signal of the external voltage or current source. Voltage voltage input Current current input

Example:

3 EPROG	
→Input	Setpoint
Si9nal	Current
L Value	0.00°C
H Value	300.00°C
Setpoint :	50.00°C

OK

OK

Example:


(i) Set the external voltage or current source output for the equivalent of a 50 °C temperature setpoint.

The value adjusted and set on the external programmer is displayed in line 4 of the LCD DIALOG-DISPLAY for inspection.

Setpoint	1:	50.00°C
IntAct	:	24.64°C
Power	:	0%
Control	:	intern

After returning the LCD display to standard display by pressing the key
 , the temperature value adjusted and set on the external voltage or current source is displayed in line 1 (Example: Eprog 50.00 °C).

This EPROG input enables the use of different voltage and current values as program parameters.

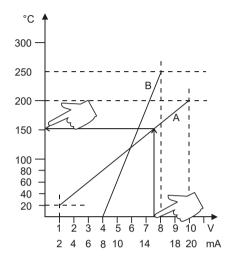

"L Value" - Setting the Low value: (See below ⊃)
1) Adjust and set the lowest desired value on the voltage or current source (Example A: 1 V).
Wait approximately 30 seconds.
2) Assign a lower temperature threshold value to this adjusted voltage/current value by pressing the appropriate keys on the keypad of the instrument (Example A: 20 °C) and set by pressing OK

"H Value" - Setting the High value: (See below ⁽²⁾)
 1) Adjust and set the highest desired value on the voltage or current source (Example A: 10 V).
 Wait approximately 30 seconds.

2) Assign an upper temperature threshold value to this adjusted voltage/current value by pressing the appropriate keys on the keypad of the instrument

(Example A: 200 °C) and set by pressing OK .

(i) Example B in the diagram illustrates that the end point values are freely selectable (e.g., 8 mA and 16 mA).

3 EPROG	
→Input	Setpoint
Si9nal	Voltage
L Value	20.00°C
H Value	000.00°C
Setpoint :	152.00°C
5ecpoint .	102.00 C
SetPoint1:	152.00°C
IntAct :	50.64°C
Power :	100 %
Control :	intern

Example out of diagram A:

.

• Adjusting the voltage source for an output of 7.6 V!

Line 5 of the LCD DIALOG-DISPLAY shows the externally set setpoint value. The instrument calculates this value from the slope of the two specified end points (in example A, 7.6 V corresponds to an external setpoint temperature of 152.0 $^{\circ}$ C).



Notice:

If this adjustment is not correctly performed at two different points, the setpoint setting will be incorrect.

Important:

The usable temperature range between > L Value < and > H Value < is limited to the configured working temperature range of the circulator or unit combination. For the working temperature range, see technical specifications.

9.9.4. External Stand-by input

2 INPUTS/0	JTPUTS
EPROG	Setpoint
→Ext.StBу	inactiv
AlarmOut	StdBy

Ext. StBy (External STAND-BY input)

Stand-By input for external switch-off.

Possible parameters:

inactive - Stand-By input is ignored

active - Stand-By input is active

Activate the standby input:

- 1. Under menu item >Ext. StBy<, set the parameter to >active<.
- 2. Connect an external contact "AK" (e.g., for external switch-off) or an alarm contact of the superordinated system.

If the connection between pin 2 and pin 3 is interrupted by opening the contact "AK", a complete shutdown of the circulating pump and heater is effected, and the unit enters the condition "**E OFF**".


As long as the contact remains open, line 4 of the LCD DIALOG-DISPLAY flashes and displays the message "Stand-By".

If the contact is reclosed, the instrument returns to the standby state and " **E OFF** " is displayed. Press **OK** to start.

(i) Additional tips for using the STAND-BY input:

The standby function can be used in conjunction with the AUTOSTART feature.

- 1. If the Autostart function is NOT turned ON, the standby input is used as described above.
- 2. If the Autostart function is enabled, the instrument will revert back to the original method of entering the setpoint (i.e., keypad, RS232, analog input, etc.).

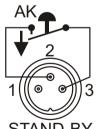

STAND-BY

Set	:	XX.XXª	С
IntAct	:	XX.XX°	C
ProgNr.	:2	Step:	2
***** 9	Stand	1By ****	*

Entering the setpoint with the keypad. e.g.,


As described above, a complete shutdown is accompanied by the "Stand-By" display and the OFF state. The programmable controller starts again when the contact is reclosed. The temperature of the bath fluid changed during the Stand-By state.

Entering the setpoint with the programmer.

The display "StandBy" appears. The setpoint value and the time are both held at the current value. The temperature of the bath fluid will be held constant at this temperature. The programmer continues once the contact is reclosed.




 Δ Notice: this is not an actual emergency switch-off.

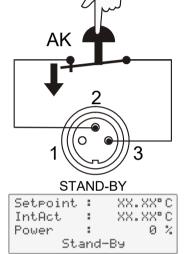



9.9.5. Alarm-output

2 INPUTS/OUT	PUTS
Ext.StBy	inactiv
→AlarmOut	StdBy

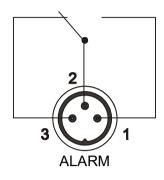

Alarm output (10)


(for external alarm signal)

This socket is a potential-free change-over contact. With the adjustments in the menu >Inputs/Outputs<, all operating conditions can be signaled without having to change the pin assignments.

3 ALARM OUTPUT	3 ALARM OUTPUT	3 ALARM OUTPUT
Function StdBy →TYPE normal	TYPE 💠 normal	TYPE 🗢 invers

Meaning of the terms under menu item >Function< :

The circulator is in condition

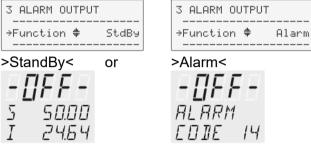

For >Type: normal<, pins 2 and 3 are connected in any case according to the selected >Function<.

For >Type: inverse<, pins 2 and 1 are connected in any case according to the selected >Function<.

Switching capacity	max.	30 W/40 VA
Switching voltage	max.	125 V~/-
Switching current	max.	1 A

10. Troubleshooting guide / Error messages





ЧП

NARNTNG

FATF

Alarm without shutdown:

Alarm with complete shutdown:

The code for the cause of alarm is indicated on the VFD COMFORT-DISPLAY. The warning signal sounds in regular intervals. The messages appear every 10 seconds.

OK to stop the signal Press the key Press OK to indicate the help text in the LCD DIALOG-DISPLAY when alarm messages appear. The circulator is operated without bath fluid, or the liquid level is insufficient. ALARM Replenish the bath tank with the bath fluid. Tube breakage has occurred (insufficient filling level due to excessive bath Low liquid level or fluid pumped out). Replace the tubing and replenish the bath tank with the float brocken. Check float and fill bath fluid. medium! The float is defect (e.g., because damaged in transit). Repair by authorized JULABO service personnel. ALARM During the self-test after switch-on a short-circuit is registered between pin 2 and pin 4 of the control cable or the control cable is interrupted during EDIE D2 operation. Control cable of the Reconnect the cable or eliminate the short-circuit. cooling machine is defective. Ckeck connection! Excess temperature warning • NABNINE AL ARM or ED 11 E - D 3 EDJE D3 Excess temperature alarm Actual temperature obove high tempera-Warn-Type: Set to ture value. >Warning< or >Alarm< Check limits Low temperature warning NABNING AL ARM or EDIE D4 EDIE D4 Low temperature alarm. Actual temperature below low tempera-Warn-Type: Set to ture value. >Warning< or >Alarm< Check limits! ALARM The cable of the working temperature sensor is disconnected or shortcircuited. EDIE OS Internal working sensor is shorted or interrupted. Call service!

Troubleshooting guide / Error messages

ALARM EDJE DE Sensor difference b.	Defect of the working or excess temperature sensor. The working temperature and excess temperature sensors report a temperature difference of more than 35 K.
working- and safety- temp. Check pump stage and viscosity!	
RLARM EDIE DI	Other errors (I ² C-BUS errors)
Internal hardware error.	
Call service! RLRRM EDIE 12	Error in A/D converter
Error in A/D converter of the measurin9 system. Call service!	
ALARM	Excess temperature sensor is defective.
EDIE IH	The safety temperature is below the working temperature setpoint. Set the
Temperature exceeds the adjusted safety- temperature. Check adjustment!	safety temperature to a higher value.
ALARM EDIE IS	External control selected, but external Pt100 sensor is not connected or is defective.
External sensor is not connected or brocken. Check the external sensor!	
WARNING EDIE 20	Cooling of the condenser is impaired. Clean the air-cooled condenser. For a water-cooled condenser, check the flow rate and temperature of the cooling water.
WARNING EDDE 21	The first-stage compressor does not work. After a short cooling interval, the compressor motor will be automatically reconnected and the message "CODE 21" will no longer appear.
WARNING	The second-stage compressor does not work.
E01E 22	<u>Cooling compressor overload protection</u> The motor of the cooling compressor is equipped with an overload protector, which will be activated in case of excessive temperature in the capsule or excessive current draw.
	Causes of motor shut-off: - poor air circulation - short distance to walls - dirt accumulated on condenser - high ambient temperature - compressor switched off and on at short intervals
WARNING EDIE 23	Excess temperature in the first-stage compressor.
WRRNING EDIE 24	Excess temperature in the second-stage compressor.

_	1151-56
WARNING EDIE 25	Short-circuit in the control cable for the cooling machine during self-test.
RLARM EDIE 33	Cable of the excess temperature sensor disconnected or short-circuited.
Internal safety- temperature sensor shorted/interrupted. Call service!	
RLARM EDIE 38	Ext. Pt100 sensor input without signal, but setpoint programming set to external Pt100.
Ext. sensor input without signal but setpoint programming set to ext. Pt100!	
WARNING EDIE 40	The early warning system for low level signals a critical fluid level. Replenish the bath tank with bath fluid.
45 ~ 4 S ~	After eliminating the malfunction, turn the mains power switch off and on again to cancel the alarm state. If the unit cannot be returned to operation, contact an authorized JULABO service center.
CONFIGURATION ERROR CONFIRM BY PRESSING : <ok></ok>	Special message "CONFIGURATION ERROR" The configuration of the circulator does not correspond to its present use. Press OK to automatically perform a one-time modification of the configuration. Then contact an authorized JULABO service center.

Disturbances that are not indicated.

The electronic pump motor is protected against overload by an electronic current limiter. If the viscosity of the bath fluid is or becomes too high, the motor stops running.



Cooling machine: Mains circuit breakers (resettable) 16 A



Fuses:

The mains fuses on the rear of the unit may easily be replaced as shown at left.

Circulator: Fine fuses (15)– T 1.25 A, 250 V~ , D 5 x 20 mm Circulator: Fine fuses (16a)– T 16 A, 250 V~ , D5 x 20 mm



Warning:

Before exchanging the fuses, turn off the mains power switch and disconnect the power plug from the mains socket!

Only use fine fuses with a nominal value as specified.

Example:

Manufacturer	Supplier	Туре	Order-No.
Schurter	Schurter	G-fuse link SPT T16A 5x20mm	No. 0001.2516
Wickmann	Wickmann	G-fuse link T1,25A 5x20 mm	No. 19195

11. Electrical connections




Notice:

Use shielded cables only.

The shield of the connecting cable is electrically connected to the plug housing.


The unit ensures safe operation if connecting cables with a maximum length of 3 m are used. The use of longer cables does not affect proper performance of the unit, however external interferences may have a negative impact on safe operation (e.g. cellular phones).

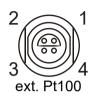

Socket for external Pt100 sensor

Pin assignment:		
<u>Pin</u>	Signal	
1	+	
2	U+	
3	U-	
4	I-	

The shield of the connecting cable is electrically connected to the plug housing and the sensor tube.

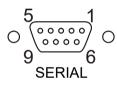

RS232/RS485 serial interface

This port can be used to connect a computer with an RS232 or RS485 cable for remote control of the circulator.

Pin assignments RS232:

Pin 2	RxD	Receive Data
Pin 3	TxD	Transmit Data
Pin 5	0 V	Signal GND
Pin 7	RTS	Request to send
Pin 8	CTS	Clear to send

Pin 1; 4; 6, 9 Reserved - do not use!

Pin assignments RS485:

В	
0 V	Signal GND
+5 V (max	. 50 mÅ)
A	
	0 V +5 V (max

Pin 1; 2; 4; 7; 9 Reserved - do not use!



Notice: RS485 serial interface

The max. load current between pin 5 and pin 6 is 50 mA.

RS232 interface cable		
Circulator (9-pol)		PC (9-pol)
Pin 2 RxD	\Leftrightarrow	Pin 3 TxD
Pin 3 TxD	\Leftrightarrow	Pin 2 RxD
Pin 5 GND	\Leftrightarrow	Pin 5 GND
Pin 7 RTS	\Leftrightarrow	Pin 8 CTS
Pin 8 CTS	\Leftrightarrow	Pin 7 RTS

Accessories:	Order No.	Description
	8 980 073	RS232 interface cable 9-pol./9-pol., 2,5 m
	8 900 110	USB interface adapter cable

✤/ Control output

The connector may only be used for control of a JULABO refrigerated circulator or JULABO MVS Solenoid valve controller for cooling water.

Pin assignment:

- Pin Signal (only with attached JULABO equipment)
- 1 +24 V (I max. current 25 mA)
- 2 0 V
- 3 Alarm relay
- 4 Reserved do not use!
- 5 Cooling pulse

(for ext


Alarm output

Stand-by input

(for external alarm signal)

This potential-free change-over contact is activated in case of an alarm when pins 2 and 3 are connected.

Switching capacity	max.	30 W / 40 VA
Switching voltage	max.	125 V~/-
Switching current	max.	1 A

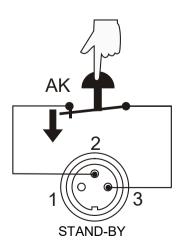


ALARM


3

(for external switch-off) Pin assignment: Pin


ssignment:	<u>Pin</u>	<u>Signal</u>
-	1	not connected
	2	5 V / DC
	3	0 V

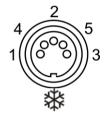

Programmer input / temperature recorder output

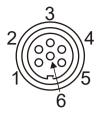


Pin		Signal		
1 Voltage output	Channel 1	0 10 V		
2 Voltage output	Channel 2	0 10 V		
3 GND for outputs	0 V			
4 Programmer input	EPROG	0 to 10 V / 0 to 20 mA		
5 Current output	Channel 3	0 to 20 mA / 4 to 20 mA		
6 GND for Progammer 0 V				

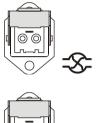

Control connectors (13, 14)

The control output (13) is not powered in the OFF condition. An external pump motor can be connected. Output voltage: 230 V~ / max. 1.25 A

The control output (14) can be adapted to solenoid valves for different tasks.

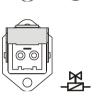

12. Remote control

12.1. Setup for remote control

$ \begin{array}{c} 5 \\ 0 \\ 9 \\ 0 \\ 0 \\ 0 \\ 0 \\ 0 \\ 0 \\ 0 \\ 0 \\ 0 \\ 0$	1.	Check the interface parameters for both interfaces (on the circulator and PC) and make sure they match. In the >Interface< menu, set the menu item >Type< to >RS232< or >RS485<.
R - FIEE -	2.	In the >Configuration< menu, set the menu item >Setpoint< to >RS232< or >RS485<.
5 5000 I 2464	3.	Connect both units with an interface cable.

Like all parameters that can be entered through the keypad, interface parameters are stored in the memory even after the circulator is turned off.

12.2. Communication with a PC or a superordinated data system

If the circulator is put into remote control mode via the configuration level, the VFD COMFORT-DISPLAY will read "R -OFF-" = REMOTE STOP. The circulator is now operated via the computer.

In general, the computer (master) sends commands to the circulator (slave). The circulator sends data (including error messages) only when the computer sends a query.

In remote control mode, the start command and all values to be set must be resent by the PC via the interface in case of a power interruption. AUTOSTART is not possible.

A transfer sequence consists of:

- address (RS485 interface only)
- command
- space (⇔; Hex: 20)
- parameter (decimal separation with a period)
- end of file (,,; Hex: 0D)
- The response (data string) after an IN command is always followed by a line feed (LF, Hex: 0A).

The commands are divided into IN or OUT commands. IN commands: retrieve parameters OUT commands: set parameters

ເສ

Important times for a command transmission:

To ensure secure data transfer, the time gap between two commands should be at least 250 ms.

The circulator automatically responds to an **in** command with a data string followed by an LF (Line Feed). The next command should be sent only after 10 ms.

(P

The **OUT** commands are valid only in remote control mode.

Examples:

When the RS485 interface is used, the three-digit instrument address precedes each command. (example: address Ad32 = A032)

Command to set the working temperature >Setpoint1< to 55.5 °C OUT_SP_00 ⇔ 55.5 J A032_OUT_SP_00 ⇔ 55.5 J

Command to retrieve the working temperature >Setpoint1< IN_SP_00با A032_IN_SP_00با Response from the circulator: 55.5با LF A032_55.5با LF



Notice: OUT commands

Temperature values can be displayed on the circulator in °C or °F. The settings sent via the interface must also be made in the selected unit according to this setting.

12.3. List of commands

OUT commands: Setting parameters or temperature values.

Command	Parameter	Response of circulator
OUT_MODE_01	0	Use working temperature >Setpoint1<
OUT_MODE_01	1	Use working temperature >Setpoint2<
OUT_MODE_01	2	Use working temperature >Setpoint3<
OUT_MODE_02	0	Selftuning "off". Temperature control using the stored parameters.
OUT_MODE_02	1	Selftuning "once" Single selftuning of the controlled system after the next start.
OUT_MODE_02	2	Selftuning "always" Continual selftuning of the controlled system whenever a new setpoint is to be reached.
OUT_MODE_03	0	Set external programmer input to voltage. Voltage 0 V 10 V

Remote control

	Response of circulator	
1	Set external programmer input to current.	
	Current 0 mA 20 mA	
0	Temperature centrel of internal beth	
	Set the control dynamics - standard	
XXX.XX	Set working temperature. "Setpoint 1"	
xxx.xx	Set working temperature. "Setpoint 2"	
XXX.XX	Set working temperature. "Setpoint 3"	
	Cot high town croture warning limit "OverTown"	
XXX.XX	-100 +100 [%]	enace
x	Set the pump pressure stage. (1 4)	
X.X	CoSpeed for external control 0 5.0.	
xxx	Xp control parameter of the internal controller.	0.1 99.9
xxx	Tn control parameter of the internal controller.	3 9999
xxx	Tv control parameter of the internal controller.	0 999
xxx	Xp control parameter of the cascade controller.	0.1 99.9
xxx	Proportional share of the cascade controller.	1 99.9
xxx	Tn control parameter of the cascade controller.	3 9999
xxx	Tv control parameter of the cascade controller.	0 999
xxx	Maximum internal temperature of the cascade co	ontroller.
xxx	Minimum internal temperature of the cascade co	ntroller.
xxx	Band limit (upper) 0 200 °C	
xxx	Band limit (lower) 0 200 °C	
-xxx	Set the desired maximum cooling power (0% to Note: Enter the value with a preceding negative This setting is meaningful only for FP cooling ma	sign!
xxx	Set the desired maximum heating power (10% to	
	0 1 0 1 0 1 0 1 0 1 ×xx.xx ×xx.xx ×xx.xx ×xx.xx ×xx.xx ×xx.xx ×xx.xx ×xx.xx ×xx.xx ×xx ×xx	Current 0 mA 20 mA 0 Temperature control of internal bath. 1 External control with Pt100 sensor. 0 Stop the unit = R –OFF 1 Start the unit. 0 Set the control dynamics - aperiodic 1 Start the unit. 0 Set the control dynamics - aperiodic 1 Set the control dynamics - standard xxx.xx Set working temperature. "Setpoint 1" xxx.xx Set working temperature. "Setpoint 2" xxx.xx Set working temperature. "Setpoint 3" xxx.xx Set high temperature warning limit "OverTemp" xxx.xx Set low temperature warning limit "SubTemp" xxx.xx Set low temperature warning limit "SubTemp" xxx.xx Set low temperature warning limit "SubTemp" xxx.xx Set the pump pressure stage. (1 4) x.x CoSpeed for external control 0 5.0. xxx Xp control parameter of the internal controller. xxx Xp control parameter of the internal controller. xxx Xp control parameter of the cascade controller. xxx Xp control parameter

IN commands: Asking for parameters or temperature values to be displayed.

Command	Parameter	Response of circulator
VERSION	None	Version number of the software (V X.xx)
STATUS	none	Status message, error message (see page 87)
IN_PV_00	none	Current bath temperature.
IN_PV_01	none	Heating power being used (%).
IN_PV_02	none	Temperature value registered by the external Pt100 sensor.
IN_PV_03	none	Temperature value registered by the safety sensor.
IN_PV_04	none	Setpoint temperature ("SafeTemp") of the excess temperature protection
IN_SP_00	none	Working temperature "Setpoint 1"
IN_SP_01	none	Working temperature "Setpoint 2"
IN_SP_02	none	Working temperature "Setpoint 3"
IN_SP_03	none	High temperature warning limit "OverTemp"
 IN_SP_04	none	Low temperature warning limit "SubTemp"
IN_SP_05	none	Setpoint temperature of the external programmer (REG+E-PROG).
IN_SP_06	none	Temperature indication in °C or °F
IN_SP_07	none	 Adjusted pump stage in the -OFF- condition. Pump stage corresponding to the effective rotational speed after start.
IN_SP_08	none	Value of a flow rate sensor connected to the E-Prog input
IN_SP_09	none	Value of a pressure sensor connected to the E-Prog input
IN_PAR_00	none	Temperature difference between working sensor and safety sensor
IN_PAR_01	none	Te - Time constant of the external bath.
IN_PAR_02	none	Si - Internal slope
IN_PAR_03	none	Ti - Time constant of the internal bath.
IN_PAR_04	none	CoSpeed - Band limit (max. difference between the temperatures in the internal bath and external system).
IN_PAR_05	none	Factor pk/ph0: Ratio of max. cooling capacity to max. heating capacity
IN_PAR_06	none	Xp control parameter of the internal controller.
IN_PAR_07	none	Tn control parameter of the internal controller.
IN_PAR_08	none	Tv control parameter of the internal controller.
IN_PAR_09	none	Xp control parameter of the cascade controller.
IN_PAR_10	none	Proportional share of the cascade controller.
IN_PAR_11	none	Tn control parameter of the cascade controller.
IN_PAR_12	none	Tv control parameter of the cascade controller.

Remote control

Command	Parameter	Response of circulator
IN_PAR_13	none	Adjusted maximum internal temperature of the cascade controller.
IN_PAR_14	none	Adjusted minimum internal temperature of the cascade controller.
IN_PAR_15	none	Band limit (upper)
IN_PAR_16	none	Band limit (lower)
IN_MODE_01	none	Selected setpoint: 0 = Setpoint 1 1 = Setpoint 2 2 = Setpoint 3
IN_MODE_02	none	Selftuning type: 0 = Selftuning "off" 1 = Selftuning "once" 2 = Selftuning "always"
IN_MODE_03	none	Type of external programmer input: 0 = Voltage 0 V to 10 V 1 = Current 0 mA to 20 mA
IN_MODE_04	none	Internal/external temperature control: 0 = Temperature control with internal sensor. 1 = Temperature control with external Pt100 sensor.
IN_MODE_05	none	Circulator in stop/start condition: 0 = stop 1 = start
IN_MODE_08	none	Adjusted control dynamics 0 = aperiodic 1 = standard
IN_HIL_00	none	Max. cooling power (%).
IN_HIL_01	none	Max. heating power (%).

12.4. Status messages

Status messages	Description
00 MANUAL STOP	Circulator in "OFF" state.
01 MANUAL START	Circulator in keypad control mode.
02 REMOTE STOP	Circulator in "r OFF" state.
03 REMOTE START	Circulator in remote control mode.

12.5. Error messages

Error messages	Description
-01 LOW LEVEL ALARM	Low liquid level alarm
-02 REFRIGERATOR ALARM	Control cable of the refrigerated circulator or MVS solenoid valve controller short-circuited or disconnected.
-03 EXCESS TEMPERATURE WARNING	High temperature warning
-04 LOW TEMPERATURE WARNING	Low temperature warning.
-05 WORKING SENSOR ALARM	Working temperature sensor short-circuited or disconnected.
-06 SENSOR DIFFERENCE ALARM	Sensor difference alarm. Working temperature and safety sensors report a temperature difference of more than 35 K.
-07 I ² C-BUS ERROR	Internal error when reading or writing the I ² C bus.
-08 INVALID COMMAND	Invalid command.
-09 COMMAND NOT ALLOWED IN CURRENT OPERATING MODE	Invalid command in current operating mode.
-10 VALUE TOO SMALL	Entered value too small.
-11 VALUE TOO LARGE	Entered value too large.
-12 TEMPERATURE MEASUREMENT ALARM	Error in A/D converter.
-13 WARNING : VALUE EXCEEDS TEMPERATURE LIMITS	Value lies outside the adjusted range for the high and low temperature warning limits. But the value is stored.
-14 EXCESS TEMPERATURE PROTECTOR ALARM	Excess temperature protector alarm
-15 EXTERNAL SENSOR ALARM	External control selected, but external Pt100 sensor not connected.
-20 WARNING: CLEAN CONDENSOR OR CHECK COOLING WATER CIRCUIT OF REFRIGERATOR	Cooling of the condenser is impaired. Clean air-cooled condenser. Check the flow rate and cooling water temperature for a water-cooled condenser.
-21 WARNING: COMPRESSOR STAGE 1 DOES NOT WORK	First-stage compressor does not work.

Installation of electronic module with analog connectors


Error messages	Description
-22 WARNING: COMPRESSOR STAGE 2 DOES NOT WORK	Second-stage compressor does not work.
-23 WARNING: HIGH TEMPERATURE ON COMPRESSOR STAGE 1	Excess temperature at first-stage compressor.
-24 WARNING: HIGH TEMPERATURE ON COMPRESSOR STAGE 2	Excess temperature at second-stage compressor.
-25 REFRIGERATOR WARNING	Error in the cooling machine.
-26 WARNING: STAND-BY PLUG IS MISSING	External standby contact is open.
-30 CONFIGURATION ERROR: CONFIRM BY PRESSING <ok> ON CIRCULATOR</ok>	The configuration of the circulator does not correspond to its present use. Press OK to automatically perform a one-time modification of the configuration.
-33 SAFETY SENSOR ALARM	Excess temperature sensor short-circuited or disconnected.
-38 EXTERNAL SENSOR SETPOINT PROGRAMMING ALARM	Ext. Pt100 sensor input without signal and setpoint programming set to external Pt100.
-40 NIVEAU LEVEL WARNUNG	Low liquid level warning in the internal reservoir.

13. Installation of electronic module with analog connectors



Caution:

Always turn off the unit and disconnect the mains cable from the power source before performing any service or maintenance procedures, or before moving the unit. Configuration, installation, maintenance, and repairs on the circulator may be performed only by qualified personnel.

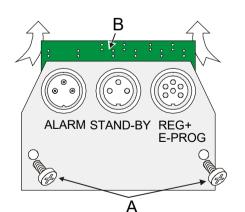

- Turn off the unit with the mains switch and disconnect the power plug.
- Remove the screws (A) and remove the dummy plate.
- Carefully insert the upper edge (B) into the rear panel from below, align the electronic module at the top and bottom, and connect the 15-pin connector by applying light pressure.
- Attach the electronic module using the screws (A).
- The circulator is ready for use. It automatically recognizes the installed electronic module.

14. JULABO Service – Online remote diagnosis

PortDef.vi
Bitte den verwendeten COM Port und die Übertragungsrate auswählen! Mit OK bestätigen!
Please choose desired COM port and the used baud rate! Confirm with OK!

COM1
4800 Baud

OK

EasyBlackBox.vi			- I X
Julabo	Ea	syBlack	kBox ersion 1.0
Einstellungen/Settings	Alarmspeicher/Alarms stored	Blackbox	
*** TEMPERATURE LIMITS Working Temperature Rang Select Temperature Range SavePoti Temperature: 90	*** : 61.00 C soint T2: 37.00 C Setpoint T3: 61.00 *** ge: -94.90 C to 200.00 C : -94.90 C to 200.00 C C : -94.90 C to 200.00 C	0C	
Speichern/Save	Hilfe/Help	Beenden/Quit	

JULABO circulators of the HighTech series are equipped with a black box. This box is implemented in the controller and records all significant data for the last 30 minutes.

In case of a failure, this data can be read out from the unit by using special software. This software is available as a free download from www.julabo.us.

Installation is easy and is performed step by step.

Please observe the instructions.


- Data read-out is possible in the conditions "OFF", "R OFF" or "ALARM".
- Connect the circulator to the computer using an interface cable.
- Start the EasyBlackBox program. The program asks for the port used (COM1,) and the baud rate of the unit. You do not have this information on hand? Simply try it out! The program continues to send the request until the correct settings are made.
- Data is read out and shown on the monitor divided into the sections
 >Einstellungen/Settings<,
 >Alarmspeicher/Alarms stored<,
 >Blackbox
 - ← see example
- After pressing >Speichern/Save<, a text file is created. The program suggests a filename ->C:\model description and barcode no.<. Modifications are possible.
- E-mail this file to service@julabo.us, JULABO's service department. JULABO is thus able to provide rapid support.

15. Cleaning / repairing the unit



Caution:

- Always turn off the unit and disconnect the mains cable from the power source before cleaning the unit.
- Prevent humidity from entering into the circulator.
- Electrical connections and any other work must be performed by qualified personnel only.

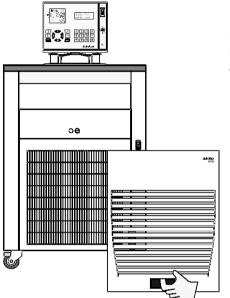

In dusty environments, the refrigerating machine's condenser should be cleaned by vacuuming every 1 to 2 months. The condenser is located behind the venting grid on the front of the unit, which can be taken off by removing the screws.

- Switch off the unit, disconnect mains power cable.
- Remove the venting grid.
- Clean the ribbed condenser with a vacuum cleaner.
- Replace the venting grid.
- Switch on the unit.

Cleaning:

For cleaning the bath tank and the immersed parts of the circulator, use low surface tension water (e.g., soap suds).

Clean the outside of the unit using a wet cloth and low surface tension water.

The circulator is designed for continuous operation under normal conditions. Periodic maintenance is not required.

The tank should be filled only with a bath fluid recommended by JULABO. To avoid contamination, it is essential to change the bath fluid from time to time.

Repairs

Before asking for a service technician or returning a JULABO instrument for repair, please contact an authorized JULABO service station.

When returning the unit:

- Clean the unit in order to avoid any harm to the service personnel.
- Attach a short fault description.
- During transport the unit has to stand upright. Mark the packing correspondingly.
- When returning a unit, take care of careful and adequate packing.
- JULABO is not responsible for damages that might occur from insufficient packing.

JULABO reserves the right to carry out technical modifications with repairs for providing improved performance of a unit.

16. WARRANTY PROVISIONS

The following Warranty Provisions shall apply to products sold in North America by Julabo ("Seller") to the entity shown as buyer ("Buyer") on Seller's invoice.

Initial Warranty

Upon Seller's receipt of payment in full for the products and subject to Buyer's compliance with the terms of sale and any other agreement with Seller relating to the products, Seller warrants to the Buyer that the products manufactured by the Seller are free from defects in material and workmanship for a period not to exceed two (2) years of operation from the date the product is shipped by Seller to Buyer (the "Initial Warranty").

EXCLUSION OF ALL OTHER EXPRESS WARRANTIES; EXCLUSION OF ALL IMPLIED WARRANTIES.

OTHER THAN THE INITIAL WARRANTY, NO OTHER EXPRESS WARRANTIES ARE MADE. ALL IMPLIED WARRANTIES OF EVERY TYPE AND KIND, INCLUDING BUT NOT LIMITED TO ANY IMPLIED WARRANTY OF MERCHANTABILITY OR FITNESS FOR A PARTICULAR PURPOSE OR USE, ARE EXCLUDED IN ALL RESPECTS AND FOR ALL PURPOSES. SELLER DISCLAIMS AND MAKES NO IMPLIED WARRANTIES WHATSOEVER.

Exclusions

The Initial Warranty does not include damage to the product resulting from accident, misuse, improper installation or operation, unauthorized or improper repair, replacement or alteration (including but not limited to repairs, replacements, or alterations made or performed by persons other than Seller's employees or authorized representatives), failure to provide (or use of improper) maintenance, unreasonable or unintended use or abuse of the product, or failure to follow written installation or operating instructions.

Buyer must return the product's record of purchase to the Seller or one of Seller's authorized representatives within thirty (30) days of the date the product is shipped by Seller to Buyer in order to make a claim under the Initial Warranty. Notwithstanding anything contained herein to the contrary, all glassware, including but not limited to reference thermometers, are expressly excluded from the Initial Warranty.

Buyer's sole remedies; Limitations on Seller's Liability

Buyer's sole and exclusive remedy under the Initial Warranty is strictly limited, in Seller's sole discretion, to either: (i) repairing defective parts; or (ii) replacing defective parts. In either case, the warranty period for the product receiving a repaired or replaced part pursuant to the terms of the Initial Warranty shall not be extended. All repairs or replacements performed by Seller pursuant to these Warranty Provisions shall be performed at one of the Seller's facility in Allentown, Pennsylvania, U.S.A. or at the facility of an authorized representative of Seller, which location shall be determined by Seller in its sole discretion; provided, however, that Seller may, in its sole discretion perform such repairs or replacements at Buyer's facility in which case Buyer shall pay Seller's travel, living and related expenses incurred by Seller in performing the repairs or replacements at Buyer's facility. As a condition precedent to Seller's obligation to repair or replace a product part under the Initial Warranty, Buyer shall (i)promptly notify Seller in writing of any such defect; (ii) shall have returned the product's record of purchase to Seller or to Seller's authorized representatives within thirty (30) days of the date the product is shipped by the seller; and (iii) assist Seller in all respects in its attempts to determine the legitimacy and basis of any claims made by or on behalf of Buyer including but not limited to providing Seller with access to the product to check operating conditions. If Buyer does not provide such written notice to Seller within the Initial Warranty period or fails to return the product's record of purchase as set forth above, Seller shall have no further liability or obligation to Buyer therefor. In no event shall Seller's liability under the Initial Warranty exceed the original purchase price of the product which is the subject of the alleged defect.

THE REMEDIES PROVIDED IN THE INITIAL WARRANTY ARE THE SOLE AND EXCLUSIVE REMEDIES AVAILABLE TO THE BUYER. NOTWITHSTANDING ANYTHING TO THE CONTRARY CONTAINED HEREIN, AND EVEN IF THE SOLE AND EXCLUSIVE REMEDIES FAIL OF THEIR ESSENTIAL PURPOSE FOR ANY REASON WHATSOEVER, IN NO EVENT SHALL SELLER BE LIABLE FOR BUYER'S MANUFACTURING COSTS, LOST PROFITS, GOODWILL, OR ANY OTHER SPECIAL, INDIRECT, PUNITIVE, INCIDENTAL OR CONSEQUENTIAL DAMAGES TO BUYER OR ANY THIRD PARTY AND ALL SUCH DAMAGES ARE HEREBY DISCLAIMED.

Assignment

Buyer shall not assign any of its rights or obligations hereunder without the prior written approval of Seller; provided, however, that if Buyer is a distributor of Seller, the rights and obligations of Buyer under these Warranty Provisions shall inure to the benefit of and be binding upon Buyer's customers who provide the product's proof of purchase to Seller pursuant to the terms set forth herein. Seller may assign any or all of its rights or obligations hereunder without Buyer's prior consent.

Governing Law

The Warranty Provisions and all questions relating to their validity, interpretation, performance, and enforcement shall be construed in accordance with, and shall be governed by, the substantive laws of the Commonwealth of Pennsylvania without regard to its principles of conflicts of law.

Waiver

Any failure of the part of Seller to insist on strict compliance with the Warranty Provisions shall no way constitute a waiver of such right. No claim or rights arising out of a breach of the Warranty Provisions by Buyer may be discharged in whole or in part by a waiver of the claim or right, unless the waiver is in writing signed by an authorized representative of Seller. Seller's waiver or acceptance of any breach by Buyer of any provisions of the Warranty Provisions shall not constitute a waiver of or an excuse for nonperformance as to any other provision of the Warranty Provisions nor as to any prior or subsequent breach of the same provision.

Freight

Seller will arrange and pay for shipping and handling for the return of the unit to the Buyer.

Out of Box Failure (OBF)

An Out of Box Failure (OBF) is defined as a product failure immediately following unpacking and installation of a newly delivered product. JULABO provides a 14-day grace period after the date of shipment, during which time the delivered product must be checked for defect. The same exclusions that apply to the regular warranty also apply to OBF classification. For example, JULABO will not be liable for transport damage, damage inflicted by the customer or any other party, or defects arising from improper installation or usage.